



Installing the bulb - BA410

30W Halogen

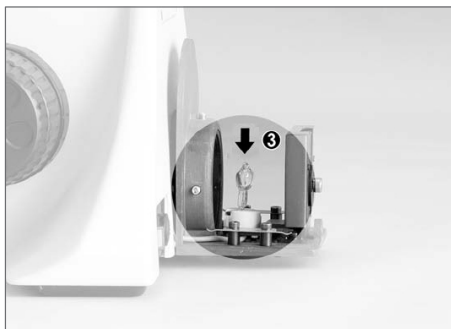


- The applicable halogen bulbs are the 6V 30W HAL high-intensity bulb.
- In order to prevent electric shock always turn the power switch off and unplug the power cord before installing or replacing the lamp.
- Loosen the lamp housing cover clamp screw using a coin and remove the cover.
- When installing the bulb, do not touch the glass surface of the bulb with bare fingers. Doing so will cause fingerprints, grease, etc., to burn onto the bulb surface, reducing the illumination provided by the bulb. If the surface is contaminated, wipe it clean using lens tissue.
- Firmly insert the bulb into the socket pinholes until it reaches the limit. Be careful not to tilt the bulb when mounting.
- Close the cover and secure it with it with lamp housing cover clamp screw.

30W LED Module

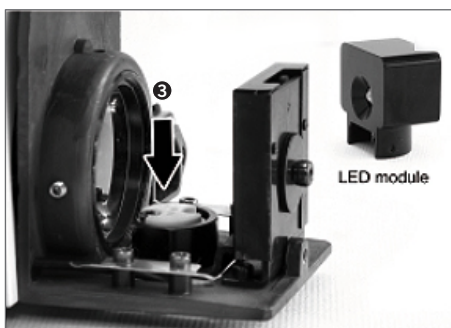

- Unplug the power supply main cable. Loosen the lamp housing cover clamp screw using a coin and remove the cover
- Firmly insert the LED module into the socket pinholes until it reaches the limit. This is a Motic patent design to exchange LED module and halogen bulb on the same socket directly.
- After the LED module installation, secure it with the clamp screw by 1.5mm hexagonal screwdriver supplied with the microscope.
- Return lamphouse cover to original position.

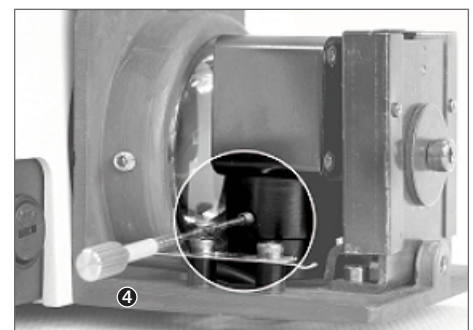



Microscope World
6122 Innovation Way
Carlsbad, CA 92009

www.MicroscopeWorld.com

PHONE: 760-438-0528

FAX: 760-438-0553

TOLL FREE USA: 800-942-0528

info@microscopeworld.com