


HG40 Parts List

Rev 2 (5/15/12)

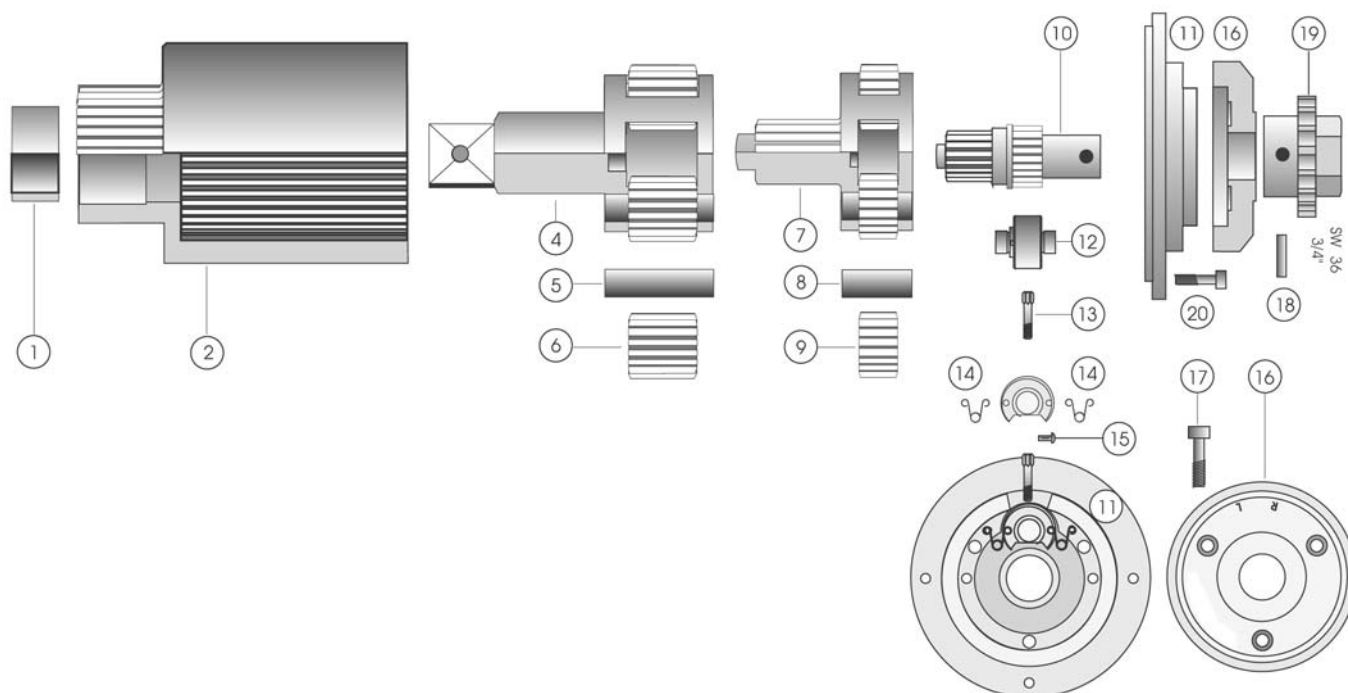

No.	Part Number	Description	Qty.
1	21-20141	Sleeve	1
2	21-10216	Gear Housing	1
4	21-10186	Gear Cage 1 1/2"	1
5	21-20014	Bearing Pins	4
6	21-10154	Planet Gears	4
7	21-10176	Gear Cage	1
8	21-20013	Bearing Pins	4
9	21-10152	Planet Gears	4
10	21-11109	Shaft	1
11	21-11103	Gear Flange	1
12	21-11111	Pawl	1
13	21-10129	Switch Screw	1
14	21-11112	Compression Springs	2
15	21-20023	Grooved Drive Stud	2
16	21-11106	Cover Plate	1
17	21-20043	Hexagon Screw	3
18	21-20022	Shear Pin	1
19	21-11114	Hand Wheel	1
20	21-20043	Hexagon Screw	4