SCHOTT ColdVision Light Source

User Manual

D20980.022 Rev D - DCN 51219

October 2019

SCHOTT North America, Inc.
122 Charlton Street, Southbridge, MA 01550-1960
Phone: (508) 765-9744, Fax: (508) 765-1299
Email: lightingimaging@us.schott.com

Website: <https://www.us.schott.com/lightingimaging/english/sensors/products/coldvision/cvls.html>

Table of Contents

Table of Contents

Safety Information	7
Introduction	7
Intended Use.....	7
Scope.....	8
Validity	8
Symbols.....	9
Related Documents.....	9
Product Related Warnings	10
Cleaning.....	11
Certifications.....	11
Assembly Location (Plant 3540).....	11
Liabilities	12
Comments.....	12
Sicherheitshinweise	13
Einführung.....	13
Bestimmungsgemäße Verwendung.....	13
Anwendungsbereich	14
Geltung.....	14
Symbole.....	15
Zugehörige Dokumente	15
Produktbezogene Warnungen	16
Reinigung	17
Zertifizierungen.....	18
Zusammenbauort (Werk 3540).....	18
Haftungen	19
Kommentare	19
Información de seguridad	20
Introducción.....	20
Uso previsto	20
Ámbito	21
Validez.....	21
Símbolos.....	22
Documentos relacionados	22
Advertencias relacionadas con los productos	23
Limpieza	24

Table of Contents

Certificaciones.....	24
Lugar de montaje (Planta 3540).....	25
Responsabilidades	25
Comentarios.....	25
Informations de sécurité.....	26
Introduction	26
Utilisation appropriée	26
Portée.....	27
Validité	27
Symboles	28
Documents connexes	28
Mises en garde relatives au produit	29
Nettoyage.....	30
Certifications	31
Site d'assemblage (Usine 3540)	31
Responsabilités	32
Commentaires.....	32
安全にお使いいただくために	33
はじめに	33
想定される使用方法.....	33
適用範囲	34
ご注意	34
シンボル	35
関連文書	35
製品に関連する警告.....	36
清掃	37
認証	37
組み立て地 (工場 3540)	37
補償	38
コメント	38
安全情報	39
引言	39
予定用途	39
範囲	40
有効性	40
标志	41

Table of Contents

相关文件	41
产品相关警告信息.....	42
清洁	43
证书	43
组装地点（第 3540 号工厂）	43
责任	44
意见	44
Light Source Operation	45
General Operation	45
Unit Operating Modes	45
Mode Selection	45
Save Current Settings.....	46
Single Channel Devices	46
Quad Channel Devices	46
Unit Mounting.....	48
Factory Resets.....	48
Soft Reset	48
Hard Reset.....	48
Reset Failures	48
Software Install	49
Windows Menu Shortcuts	50
Driver Install.....	50
CV-LS Dashboard.....	51
Windows Menu.....	53
Connections	54
Finder Requirements	55
System Status.....	56
Configurations (ADMIN Level)	58
Controls (USER Level).....	59
LED Channel Settings	60
Continuous Strobe	62
Triggered Strobe	64
Equalizer.....	66
Fan Control.....	68
Upgrade (ADMIN Level)	69
Settings (ADMIN Level)	70
General.....	71

Table of Contents

Network	72
Sockets	74
UART.....	75
Users (ADMIN Level)	76
Add / Edit User	77
Delete User	77
Diagnostics (ADMIN Level).....	78
Sockets	78
Legacy Socket.....	79
Socket Connection	79
Command Structure.....	80
Negative Acknowledgements	80
Commands	80
Binary Socket.....	81
Socket Connection	81
Fletcher16 Checksums	81
Encoding a Packet	82
Decoding a Packet.....	82
Commands	82
Product Specification	83
Environment	83
Operational Temperature	83
Storage Temperature.....	83
Humidity.....	83
Electrical.....	83
Input.....	83
Redundant Power Inputs	83
Mechanical.....	84
Weight (Typical)	84
Physical Dimensions (Inches)	84
Front Panel Description	85
Rear Panel Description.....	86
Multiport Pin Details.....	87
LED Information	88
A20980/6000K	88
A20980/3000K	89
A20980/RGBW	89
Replacement Parts.....	90

Table of Contents

Accessories.....	90
Return Material Authorization (RMA)	91
Troubleshooting.....	91
Technical Support	91
Service and RMA Policy.....	92
Warranty and Liabilities	92
RMA Request Form	93
Customer Support.....	94
Downloads	94
Technical Support	94

Safety Information

Introduction

Read these instructions carefully and look at the equipment to become familiar with the device before operation. The following special messages may appear throughout this manual to warn of potential hazards:

CAUTION

CAUTION indicates a potentially hazardous situation that, if not avoided, **can result** in injury or equipment damage.

Electrical equipment should only be installed, operated, serviced, and maintained only by qualified personnel. SCHOTT North America assumes no responsibility for any consequences arising out of the use of this material.

Intended Use

The ColdVision Light Source (CV-LS) is a high intensity LED light source intended for use with fiber optic illuminators. This light source has state of the art light output and a simplified user interface with versatile remote control abilities. The CV-LS is compatible with all standard ColdVision Series light guides. Optimal performance is achieved when coupled with larger bundles diameters, 9-14mm.

The CV-LS currently has the following options available:

- A20980/6000K (Cool White Light Source)
- A20980/3000K (Warm White Light Source)

Additional configurations available (minimum quantities and lead times may apply):

- A20980/RGBW (RGBW Light Source)
- A20980/400NM (400NM UV Light Source)

All of these product models have industry leading light outputs, with the Cool White Light Source model topping 150W EJA halogen lamp levels. Solutions with other wavelengths, including four-channel combinations, can be produced to match specific customer application requirements.

The CV-LS has many advanced features including:

- Ethernet, RS232, and USB (Virtual RS232) [Backwards compatible with the LLS 3]
- Ethernet [Robust binary interface for accelerated communications]
- Analog Controls [0-5V standard input]
- Digital Controls [0-24V standard input]
- Triggered or Continuous Strobe [up to 10KHz]
- 25 microsecond rise time on strobe pulses
- Auto sensing inputs with control lockouts via command interfaces
- Web interface for control, user management, remote upgrade, etc.

Scope

This manual is designed to provide users with the technical information needed to operate the CV-LS. The units described in this manual are:

- A20980/6000K (Cool White Light Source)
- A20980/3000K (Warm White Light Source)
- A20980/RGBW (RGBW Light Source)

Validity

SCHOTT North America cannot be held responsible for any errors in this manual. Please contact us if you have any suggestions for improvements, modification, or corrections.


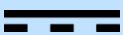
No part of this document may be reproduced in any form, or by any means, without prior authorization of SCHOTT North America.

The data, photos, and illustrations within this document do not in any way imply a contractual obligation of any kind. SCHOTT North America reserves the right to modify our products in line with our policy of continuous development. The information in this document may be modified without notice and should not be interpreted as binding on SCHOTT North America.

Safety Information

Symbols

Symbols used on this product or in this manual are described in the following table:

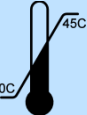Symbols	Description
	Read the manual before operating the unit.
	Caution: Surface may be hot.
	Certified by Intertek. See Certification section on page 11 for compliance information.
	European Declaration of Conformity
	RoHS Compliant, also known as Directive 2011/65/EU
	Compliant with the Waste Electrical and Electronics Equipment/Restriction of the Use of Certain Hazardous Substances in Electrical and Electronics Equipment (WEEE/RoHS) Recycling directives
	Operational Temperature Range: 0°C – 45°C
	Operational Humidity Range: 5% - 95% Noncondensing
	DC Power Input

Related Documents

Title of Document	Reference Number
CV-LS Data Sheet	
CV-LS Programming Guide	

Product Related Warnings

All relevant local safety regulations must be followed when installing and using this product. For safety reasons, and to ensure conformity with assembly standards, only the manufacturer is authorized to repair the components.

CAUTION

Never look directly at the LED output when on. LEDs have a very high luminance that can cause eye injury. Do not look at beam with optical instruments.

Use Ultraviolet (UV) protective goggles when using a UV light source.

CAUTION

Light absorbing materials have the physical property of converting incident light into heat. Damage may occur to heat sensitive or flammable light absorbing materials. To avoid such thermal damage and the potential danger of fire or burns, please adhere to the following instructions:

- Do not cover the light guide exit.
- Do not place any part of your body in the light output path.
- When illuminating heat sensitive or flammable light absorbing objects (e.g. in microscopy), special care must be taken to keep an appropriate distance between the light guide and the object(s). Special care must also be taken to select a suitable lamp brightness level in order to avoid thermal damage to the object(s).
- To avoid unnecessary strain on biological tissue caused by illumination with visible light, reduce the brightness and duration of illumination to the absolute minimum level required.
- Do not cover the front or rear vents when operating the unit. Airflow must be unrestricted to allow full light output and maintain life of the LED.
- Use only power cords with appropriate connectors for secure use.
- Disconnect power cord when unit will be left unused for long periods.
- Do not use this unit near any liquids, or in an area with excessive moisture.
- Do not place flammable materials on or near the unit.
- Use only standard glass cleaners. Do not use solvents, petroleum distillates, or any volatile or flammable liquids or foams.
- Never spill liquid on the unit.
- Do not service the unit beyond what is described in this manual. Return the light source to an authorized service center. For a Return Material Authorization (RMA), see page 91.

Cleaning

Both the housing and the front and back plates have a durable finish that should retain its original luster for many years. Cleaning the exposed areas with a commercial glass cleaner will help to maintain the finish.

Before beginning, remove the cord from the light source. Wipe the exposed areas of the housing with either a soft cloth or a paper towel moistened with a commercial glass cleaner.

CAUTION

Do not use detergents, excess water, treated cloths, harsh cleaning or flammable agents or sprays. If fluid spills into the interior, let the unit dry thoroughly before using. Periodically, any visible dust accumulation should be removed using a vacuum or commercially available can of compressed air.

Certifications

The ColdVision Light Source conforms to the provisions of the following European directives:

- 2006/95/EC (Low Voltage Directive)
- 2004/108/EC (EMC Directive)
- 2011/65/EU (RoHS)

The illumination systems observance of the basic requirements set forth in the directives mentioned above is verified by technical documentation and in full compliance with the standards listed below:

Category	Standards
EMC	IEC 61326-1:2012 FCC CFR47 Part 15B:2015 Japan Deviations
Safety	CENELEC EN 61010-1:2010 UL 61010-1:2012 CAN/CSAC22.2#61010-1:2004
Photo-biological	IEC 62471:2006

Assembly Location (Plant 3540)

SCHOTT de México, S.A. de C.V.

Carretera México-Veracruz Km.349 Venta Parada,

Amatlán de los Reyes, Veracruz México 94946

Phone: +52 (271) 7166-316

Website: www.schott.com/pharmaceutical_packaging

Safety Information

SCHOTT
glass made of ideas

Liabilities

Any warranty implied under State Law shall be limited to one year from original date of delivery to the original purchaser. Specifically excluded from SCHOTT liability is damage resulting from acts of any deity, malicious mischief, vandalism, riots, wars, improper installation, or neglect in the operation or maintenance of the unit or misunderstanding of the properties of the unit. Under no circumstances shall SCHOTT be liable for consequential or other damages of any kind, losses or expenses in connection with, by reason of the use, or inability to use, of this unit for any. The stated warranty provides the purchaser with specific legal rights, which vary from state to state. Some states, for example, do not allow exclusion of consequential damage.

Comments

We welcome your feature requests and comments about this document.

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Light Source Operation

General Operation

The CV-LS has a simplified user interface for easy startup and operation. To use the light source in normal operation, press the Power Switch and adjust the intensity control knob.

The Digital Input pins on the rear Multiport can be used to strobe the LED in 200 μ s increments. This pin can be controlled with just a contact closure switch between the Digital Input pin and GND pin, or driven low from a digital control pin. The default behavior of this pin is Active Low. When the pin is driven low or connected to GND, the LED will shut down.

More advanced controls over the light source are available through any of the digital interfaces, and are documented in the Technical User Guide (See the Downloads section on page 94).

CAUTION

Because of eye hazards, units outside the visible spectrum require a contact closure on the Digital Input pins of the Multiport to turn the LED on. The user is required to supply sufficient interlocks on this pin to protect the user from exposure to UV and Infrared (IR) light.

Unit Operating Modes

Mode Selection

There are operational modes that can be easily selected by only using the front knob and power button on the unit. These modes allow a simplified use of the unit without the more advanced techniques described in the Technical User Guide (See the Downloads section on page 94). The button on the front of the unit, when held down, has the following sequence:

- 0 to 2 seconds: No Effect (LED ring light will remain in the same state)
- 2 to 6 seconds: Mode Select (LED ring light will flash ten times per second)
- 6 to 10 seconds: Save Current Settings (LED ring light will flash two times per second)
- >10 seconds: Soft Reset (LED ring light will flash one time per second)

The mode selections are different depending on the type of device you have. See the other sections on Single Channel Devices and Quad Channel Devices below.

Light Source Operation

Save Current Settings

Once you have the state of the device to a point you would like to restore on power up, hold the button for 6 to 10 seconds to “Save Current Settings”. Upon next power up of the unit, it will return to the last saved settings.

Single Channel Devices

Single channel devices, such as the Cool White (A20980/6000K) and Warm White (A20980/3000K), have three operational modes:

- **Home (0% knob position):** This mode limits the fan to its minimum speed, and the LED output to 70% of maximum. This is the default mode, and can be used in any setting where low noise and vibration from the fan is critical. In this mode, the unit should not be used in an ambient temperature above 25°C. Operating in this mode only, the unit complies with Class B Radiated Emissions limits.
- **Industrial (50% knob position):** This mode limits the fan to 25% of its maximum speed, but increases light output to 100% of maximum. This mode can be used when more light output is needed, but with slightly higher fan noise and vibration. In this mode, the unit should not be used in an ambient temperature above 25°C.
- **Heavy Industrial (100% knob position):** This mode allows the fan to run at 100% of its maximum speed, and allows light output to 100% of maximum. This mode offers the most cooling of the LED, and will allow the unit to operate up to its maximum ambient operational temperature of 45°C. It will also extend the life of the LED at cooler ambient temperatures.

Quad Channel Devices

Quad channel devices, such as the RGBW (A20980/RGBW) have six operational modes. The mode descriptions below will be referencing the RGBW units, but can apply to any custom combination of four channels. Channel 1 is selected by default.

- **Common Channel (0% knob position):** This mode sets a scalar that is applied to all four channels. This mode is used to dim a channel mixture set by the other modes.
- **Channel 1 (20% knob position):** This mode controls the Red channel and sets its intensity.
- **Channel 2 (40% knob position):** This mode controls the Green channel and sets its intensity.
- **Channel 3 (60% knob position):** This mode controls the Blue channel and sets its intensity.
- **Channel 4 (80% knob position):** This mode controls the White channel and sets its intensity.
- **Demo (100% knob position):** This mode cycles through every combination of Channels 1-3 (Red, Green, and Blue) to demonstrate all the possible colors that the unit can generate.

Light Source Operation

Single-Color Control Example

One common use of the RGBW unit is a single color light source, where the user only needs a single color at a time, but may need to change between colors. Use the following sequence to achieve this type of operation:

1. Press and hold the button for more than ten seconds. The LED ring light will be blinking once per second.
2. Release the button to return to factory defaults.
3. Default settings will put the common channel at 100% and the knob will control channel 1 (Red).
4. Press and hold the button for two seconds. The LED ring light will start blinking ten times per second.
5. Once the LED ring light is blinking ten times per second, turn the knob to the desired channel. See the available modes above, IE 40% position for Green.
6. Release the button before six seconds has elapsed. This will change the channel and the corresponding color will turn on. **Note: If the button was held for more than six seconds, the LED ring light will start blinking two times per second. Releasing the button at this time will only save the current settings, not change the channel.**
7. At this point, the unit will act as a single channel light source of the selected color.
8. You can press and hold the button for six seconds (the LED ring light will be flashing two times per second), then release to store this operating mode.
9. To change to a different color unit, turn the knob to 0% to disable the current channel.
10. Follow steps 4 through 7 to select a new channel.

Multi-Color Control Example

A second common use would be to set a color mix using multiple channels, then dimming the pre-mixed color. Use the following sequence to achieve this type of operation:

1. Follow the instructions to set up the unit for a single color.
2. Set the knob at the intensity for that color to mix with another.
3. Follow steps 4 through 7 of the Single-Color Control Example section above to select a new channel (the first channel should still be active at the intensity you last set).
4. Repeat for as many channels as you would like to mix.
5. Once the target color mix is achieved, set the unit back to the common channel (0% point of the knob) to dim the color with the knob.

Unit Mounting

The CV-LS comes pre-installed with four rubber feet on threaded posts. These are suitable for using the unit on a flat surface to provide vibration reduction and keep the unit from slipping. These feet are secured using lightweight thread-locker to keep them from unscrewing during use.

To mount the unit using these threaded holes, the feet can be removed without damaging the unit. It is normal, due to the thread-locker, for removal of the feet to require some amount of force.

Factory Resets

There are two forms of factory resets available. Both can be accessed through different methods.

Soft Reset

The soft reset will restore factory settings to the unit. This can be achieved by first applying power to the unit, then holding the Power Switch for 10 seconds. Upon release, the unit will restore factory settings and reboot. Note: This will also reset network settings.

The soft reset can also be accessed through the digital interfaces and is documented in the Technical User Guide (See the Downloads section on page 94).

Hard Reset

The hard reset will restore factory firmware and factory settings. This can be achieved by first removing power to the unit. While holding the Power Switch, apply power to the unit, and continue holding the Power Switch for five seconds.

The Indicator light will come on for a five-second interval. Release the Power Switch at any time within this interval for the unit to perform a hard reset.

If the Indicator starts flashing twice per second, the five-second interval has elapsed. Go back to the beginning of the Hard Reset section.

Once the Indicator light starts flashing 10 times per second, the unit is restoring the factory firmware. When the Indicator light stops flashing, the hard reset is complete.

Reset Failures

If at any point you have issues restoring the unit to factory settings, see the Return Material Authorization (RMA) section on page 91 for additional instructions.

Software Install

Software Install

Download the full installer package from the Downloads Page (See the Downloads section on page 94).

This install includes:

- CV-LS User Manual (this document)
- CV-LS Data Sheet
- CV-LS Windows USB Driver
- CV-LS 3D Models
- CV-LS Firmware
- CV-LS Dashboard
- CV-LS Dashboard Source Code
- CV-LS Programming Guide

The default install location for these items is:

C:\SCHOTT North America, Inc\CV-LS, ColdVision Light Source\

Note: This location can be changed during the install, but should not be placed in a **System Folder**. Placing these files in a **System Folder** will prevent the application from accessing its settings files.

Windows Menu Shortcuts

Once the package is installed, you will find shortcuts in your start menu, as shown in Figure 1 below.

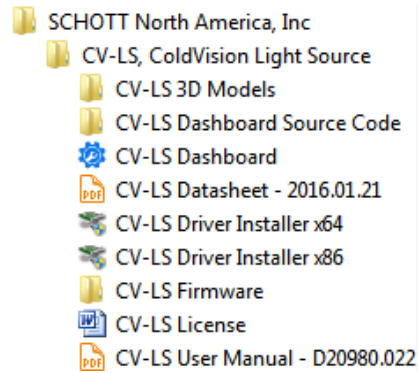
Figure 1: Windows Start Menu Shortcuts

File/Folder Name	Description
CV-LS 3D Models	This folder contains IGES, STEP, and SolidWorks Part Files.
CV-LS Dashboard Source Code	This folder contains the full C# .net source code used to create the CV-LS Dashboard Application. This code is released under the MIT Open Source License. This provides source code the user can reference, or use parts of in their own applications, to quickly interface with the CV-LS units.
CV-LS Dashboard	CV-LS Control Application. This application can fully control the CV-LS units over USB, UART, or Ethernet.
CV-LS Datasheet	Datasheet for the CV-LS product line.
CV-LS Driver Installer (x64, x86)	These applications will install the USB Virtual RS232 drivers into windows. Administrative Rights Required.
CV-LS Firmware	This folder contains the firmware file used to upgrade units through the web interface.
CV-LS License	The CV-LS Dashboard is released under the MIT Open Source License - the text of which is contained in this file.
CV-LS User Manual	User Manual for the ColdVision Light Source.

Driver Install

Install the SCHOTT CV-LS Install Package (See the Software Install section on page 49). The driver installers can be accessed from the start menu links, or directly from the install folder. The default location is:

C:\SCHOTT North America, Inc\CV-LS, ColdVision Light Source\CV-LS Driver

This folder contains a simple installer for x86 and x64 architectures, as well as the signed driver files for manual installs. **Administrative Rights** are required to install the drivers.

CV-LS Dashboard

The CV-LS Dashboard Application has many features that can be used to control any CV-LS. Some advanced settings are currently available only through the web interface, which can be launched from the Dashboard.

There are many features built into the Dashboard that are beyond the scope of this manual. Most are self-explanatory, while some may be covered in more detail in Webinars.

Each section of the Dashboard corresponds to a section of the web interface. In the following sections, there will be a screenshot from both the Dashboard and Web Interface. Descriptions of features will indicate from which interfaces they are accessible.

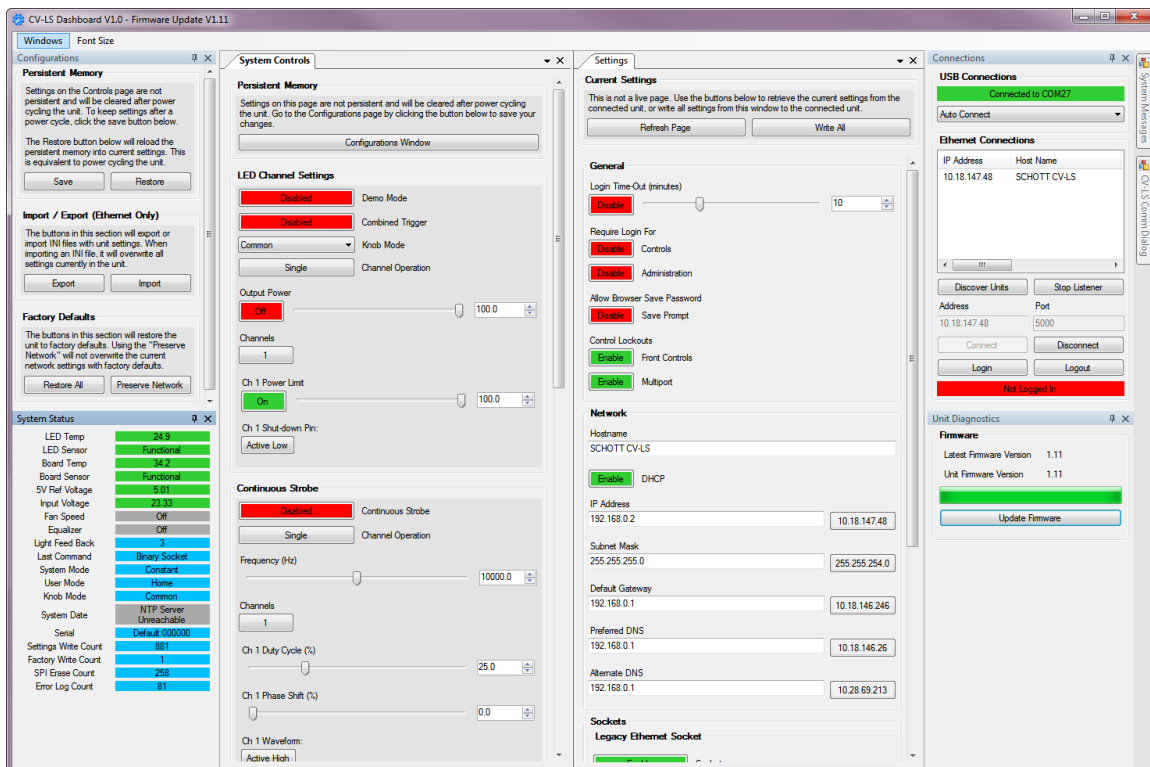
Figure 2: CV-LS Dashboard

All windows in the Dashboard are dock-able. If a window is closed, it can be reopened from the “Windows” menu. Windows can also be undocked from the dashboard into a separate window. Multiple undocked windows can be docked in the same external window. This can be seen in Figure 3 below.

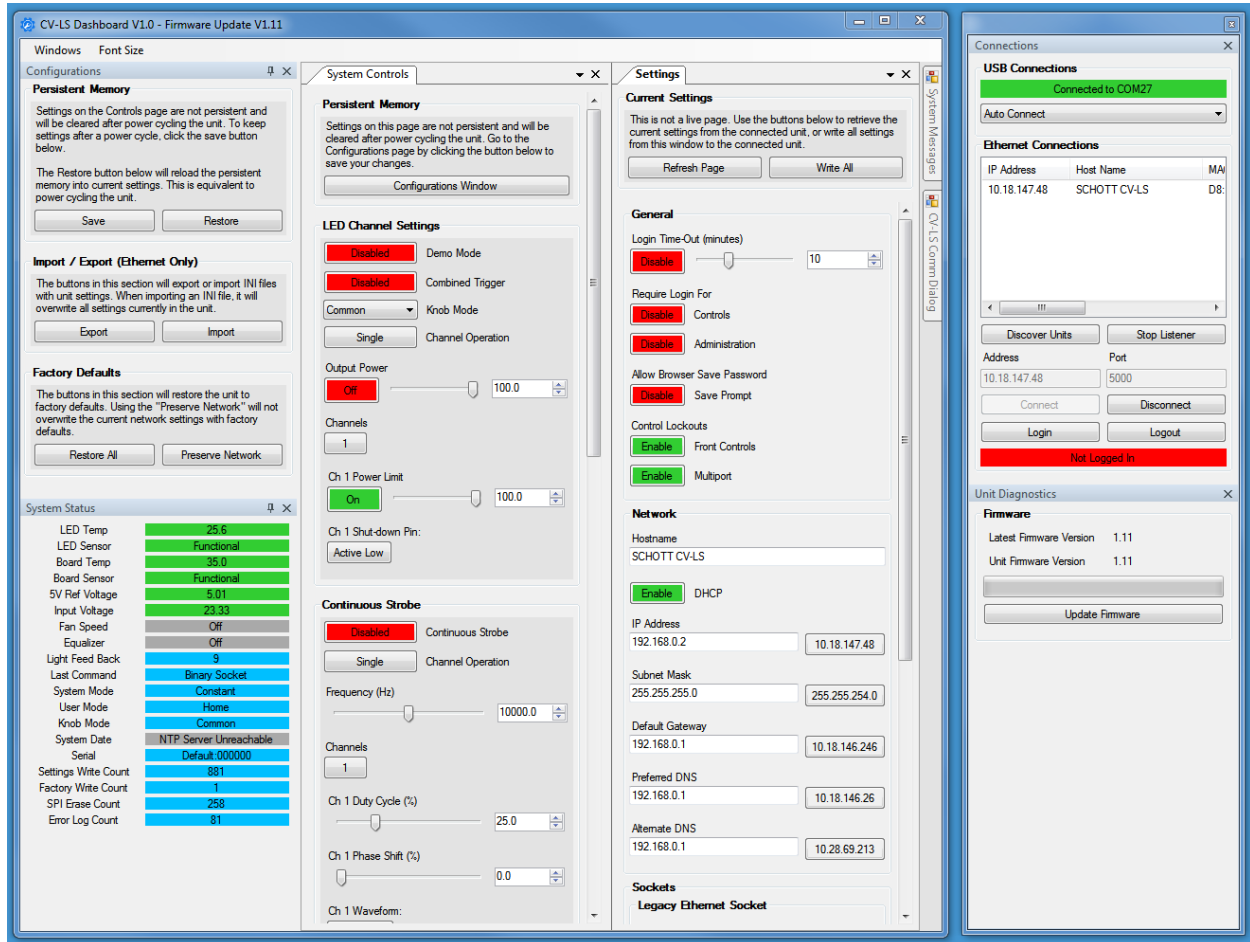
Figure 3: Controls Docked in External Window

Layouts can be saved and restored with options under the “Windows” menu, as well as reset to the default layout. Most window functions are available to all connection types, but some of the more advanced functions are available only over Ethernet connections. Advanced functions include all graphing windows, live system status, firmware upgrade, and INI import/export.

Windows Menu

The Windows Menu allows the user to open windows that are currently closed. This menu also allows the user to save / load / reset the interface.

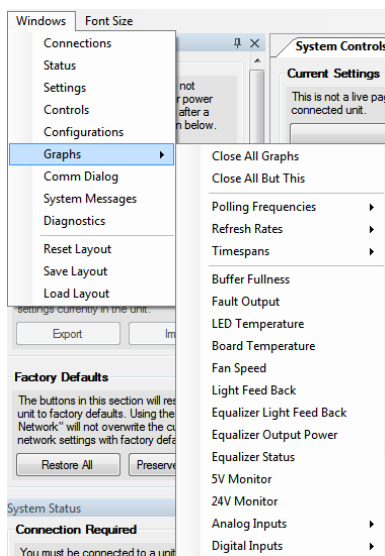
Figure 4: Window Menu

Graphing functions are highly dependent on the speed of the computer being used. The default polling rate of the graphs is 100Hz. This can be increased to 5kHz, but is not recommended unless either a very powerful computer is being used or only a few graphs are being displayed. If the graphs become backlogged with data, the connection will reset.

Note: The graph options under the Windows menu will affect all graphs in the application. You can change the settings of individual graphs by right clicking in the graph.

Color	Description
Connections	Displays the connection options and status for both USB and Ethernet.
Status	Shows the live status of the system.
Settings	ADMIN level login. Contains settings for: Login, Controls Lockout, Network, Socket, and UART.
Controls	USER level login. All LED controls for the system.
Configurations	ADMIN level login. Memory management functions.
Graphs	Contains all graphing windows. Ethernet connections only.
Comm Dialog	Telnet terminal to enter commands manually. For USB connections only.
System Messages	Any output from the Dashboard application the user may need to see for diagnostic purposes.
Diagnostics	Unit firmware upgrade.
Font Size	Allows the user to change the font size of the Dashboard Application.

Connections

The Connections section, which is specific to the Dashboard Application, is shown below in Figure 5. This section is used to connect the Dashboard to units over USB, UART, or Ethernet. If a unit is found on the network, its IP Address can be double-clicked to open the Web Interface for that unit.

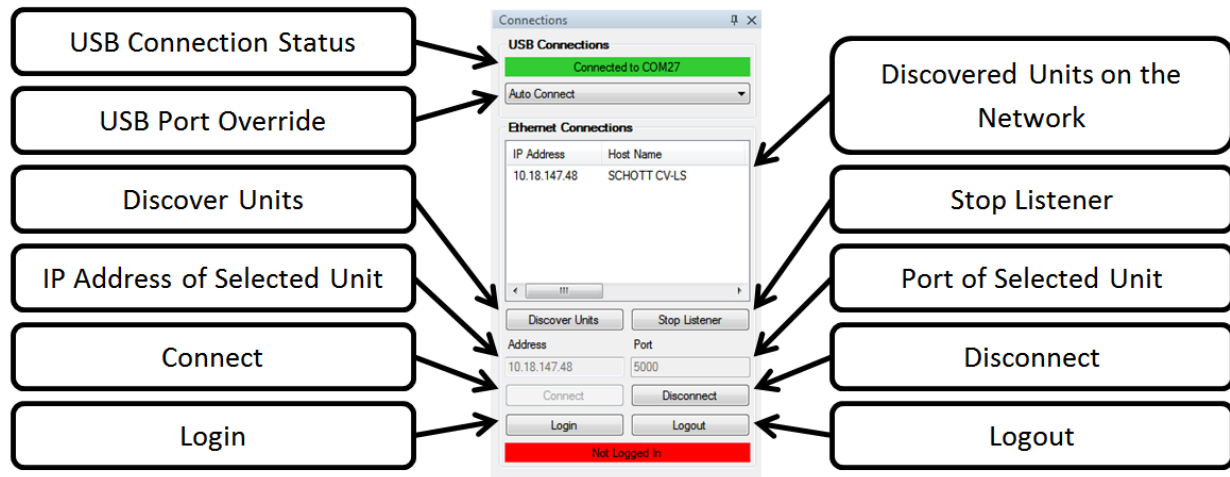
Figure 5: Connections

The Dashboard will automatically connect to the first unit plugged into USB. If a second unit is plugged into the computer, control can be switched to the second unit by changing the “USB Port Override” from “Auto Connect” to the port number of that unit. This method can also be used to connect to a unit over RS232.

The Ethernet Connections section uses a broadcast UDP packet on port 30303. Units on the network may only be listened for by one application on a computer at a time. To connect to multiple units on the network in separate applications, click “Stop Listener” in one application after connecting to a unit. Then, click “Discover Units” in the next window.

Finder Requirements

It is important to note that this method will only work if the following are true:

- 1) Broadcast UDP packets are allowed
 - a) Some networks do not allow this. Check with your IT department to see if this is blocked.
- 2) Port 30303 is not blocked
 - a) Firewalls will block the finder application. You will need either to add an exception to the firewall, or temporarily disable the firewall to discover a unit.
- 3) Unit and computer have compatible IP addresses.
 - a) The default IP settings of the unit are:
 - i) IP: 192.168.0.2
 - ii) Subnet Mask: 255.255.255.0
 - iii) Default Gateway: 192.168.0.1
 - iv) DHCP: Enabled
 - b) Both the computer and the unit should be on the same network with a DHCP server, or the computer should be set to a static IP like 192.168.0.100.

If all of these steps are followed, and no units are discovered on the network, a USB connection to the unit will display its network settings, which can then be used to diagnose connection issues.

Note: A direct Ethernet connection to a known unit IP address is possible, even if the finder application cannot find the unit on the network.

Once units are found on the network, their IP addresses will be displayed in the Ethernet Connections box. Clicking on the line once will copy the IP address into the “IP Address of Selected Unit” box. Once this is complete, clicking the “Connect” button will connect to the unit. Double-clicking on a line will open the web interface for that unit.

Certain functions of the interface can be locked out, requiring a USER or ADMIN level login. When connecting to a unit by Ethernet, these sections will be hidden with a message that they require a login of a certain level. You can login by clicking the “Login” button. At any time, you can click “Login” again to change the account that is logged in. You can also click “Logout” to logout of the current account. The bottom box will show which account is currently logged in.

CV-LS Dashboard

System Status

The System Status section shows live information about the unit. Item status is indicated by the following colors:

Color	Description
Green	Normal Operational Status.
Yellow	Warning range. Unit is still operating normally, but is near the Error threshold.
Red	Error Status. LED will not turn on if any errors exist. Fault output line will pull low (GND) if any errors exist.
Blue	Information Status.
Grey	Off status.

Figure 6: System Status (Left: Dashboard, Right: Web Interface)

Control Name	Description
LED Temp	Temperature of the LED. Normal Operational Range: 0 – 80°C Warning Range: 80 – 90°C Thermal Shutdown: 90°C
LED Status	Green when temperature sensor is fully operational. Yellow when temperature is still accurate, but sensor is degraded. Red when temperature is no longer accurate.
Board Temp	Temperature of the LED. Normal Operational Range: 0 – 65°C Warning Range: 65 – 80°C Thermal Shutdown: 80°C
Ref Voltage	5V output voltage measurement.
Input Voltage	Measurement of the current input voltage. Unit will run on 18-28V.
Fan Speed	RPM of the unit fan. Range from 7 – 23K
Equalizer	Status of the equalizer when enabled.
Light Feedback	Current light output value. Value has no units and is not calibrated. Used internally to stabilize output through equalizer functions.
Last Command	Shows the last interface that received input.
System Mode	Currently active unit mode.
User Mode	Last user selected mode.
Knob Mode	Last user selected knob mode.
System Date	When unit has internet access, time will be set from NTP servers.
Serial	Unit model and serial number.
Settings Write Count	Number of times the user has stored settings in the unit. Max value >40e6
Factory Write Count	Number of times the factory has written default settings. Max value >1.6e6
SPI Erase Count	Number of times the unit firmware has been written. Max value > 1e4
Error Log Count	Number of times a system exception has occurred. Max value > 25.6e6

Configurations (ADMIN Level)

The Configurations section is used to modify the unit memory. Current settings can be saved / restored, or reset to factory defaults in this section. All settings can also be stored as a standard format INI file, and imported into other units. These features can be accessed from both the Dashboard, as well as the Web Interface.

Note: Keep in mind that settings changed in the Controls section (page 58) are not persistent until saved in the Configurations section. This allows the user to test new settings without losing those previously saved. Once the settings are saved, the unit will return to those settings upon powering up.

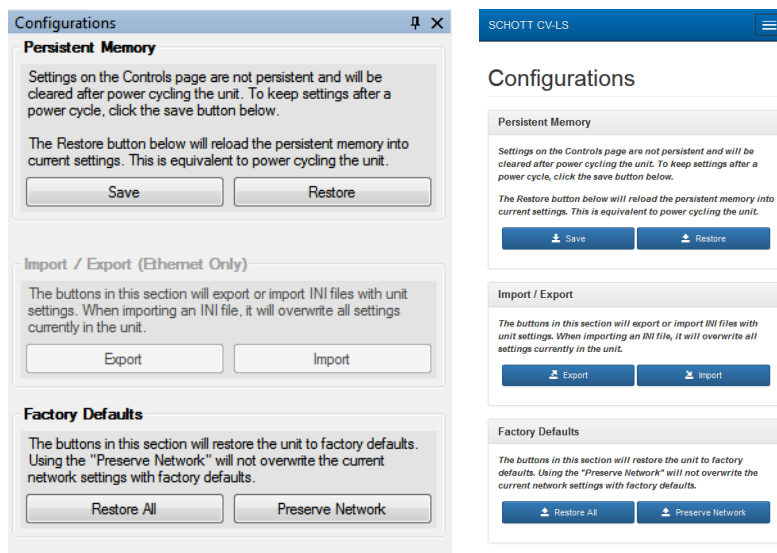
Figure 7: Configurations (Left: Dashboard, Right: Web Interface)

Control Name	Description
Save	Stores the current system settings to memory.
Restore	Restores the last saved settings to the system.
Export	Exports the current settings to an INI file.
Import	Imports an INI file into the current system settings.
Restore All	Restores all factory settings to current memory.
Preserve Network	Restores all factory settings, except network settings, to current memory.

Controls (USER Level)

The Controls section contains all of the settings pertaining to how the LED operates. There are a few differences in how the Dashboard and Web Interface behave. When the Dashboard is only connected via USB or RS232, the Controls page is static. To see the settings currently in the unit, the user must click the “Refresh Page” button at the top of the window.

Note: The “Refresh Page” button will only be shown when the Dashboard is connected to the unit by USB only.

If the Dashboard is connected via Ethernet, or if using the Web Interface, the window is continuously updating itself. For instance, if a user were to change a setting from the Web Interface, those changes would be immediately reflected in the Dashboard.

Note: There is also a message at the top of both the Dashboard and Web Interface, reminding the user that changes on these controls are not “Persistent” - meaning they will not be kept after a power cycle. To keep the changed settings, the user must manually save them from the configurations page.

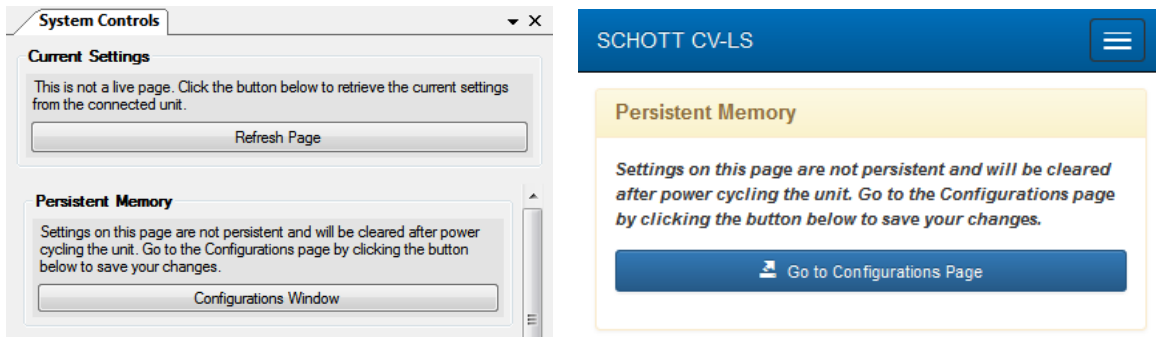
Figure 8: Controls (Left: Dashboard, Right: Web Interface)

All of the advanced operating modes for the LED can be accessed from this section. They are visually broken down into sub-sections within the interfaces. Each sub-section will be explained below.

Note: The CV-LS is a four-channel device. Each channel can be controlled independently of the others. Some of the features in the table below will not be needed in a single channel device.

LED Channel Settings

The LED Channel Settings section controls the basic functionality of the LED. When operating in this mode, the unit behaves as a simple analog dimming light source.

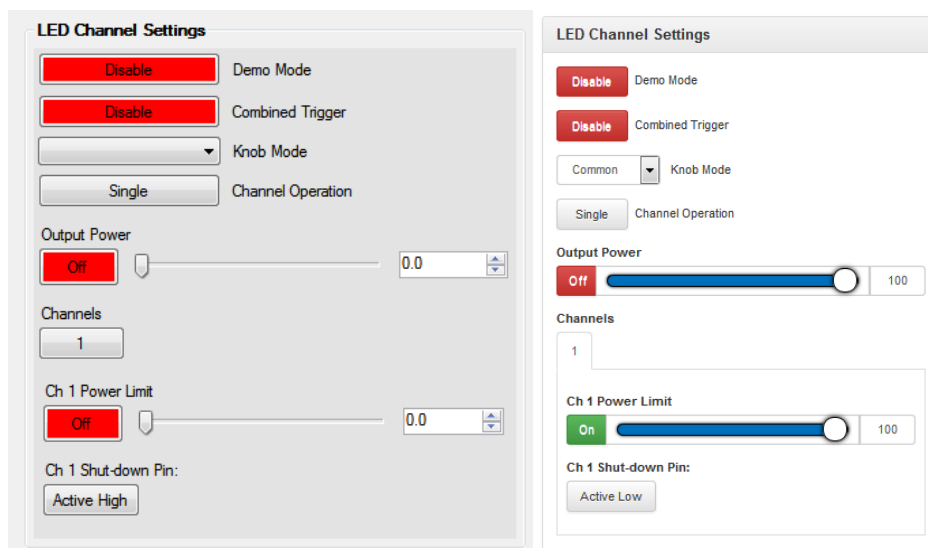
Figure 9: LED Channel Settings (Left: Dashboard, Right: Web Interface)

Control Name	Description
Demo Mode	Options: Enabled, Disabled When Demo Mode is enabled, channels 1-3 will cycle through all combinations of intensities. This function is only used for demonstrating the color mixing capabilities of the RGBW model.
Combined Trigger	Options: Enabled, Disabled When enabled, a digital input into any channel port is equivalent to applying the same digital signal to all four ports.
Knob Mode	Options: Common, Channel 1-4, Demo The front knob of the unit will control either the Common Output Power slider, or the Power Limit slider of the selected channel. Generally used to control the RGBW unit as a single color light source.
Channel Operation	Options: Single, Quad When Single is selected, the settings of Channel 1 are applied to all four channels. When Quad is selected, each channel is controlled by the respective settings for that channel.

Control Name	Description
Output Power	<p>Options: On, Off</p> <p>This button controls the LED. This must be set to ON before the LED can turn on. When control is set to ON, the LED can still be off by using the shutdown pin on the Multiport connector.</p> <p>Note: This setting is also controlled by the button on the front panel of the unit.</p> <p>Note: If this setting is ON, but the LED is not coming on, check the individual channel settings.</p>
Output Power Slider and Numeric Up Down	<p>Range: 0-100.0</p> <p>This sets the output power of the LED.</p> <p>Note: This setting is also controlled by the knob on the front panel of the unit.</p>
Channels	<p>Options: 1, 2, 3, 4, All</p> <p>When in Quad mode, all options are displayed to allow the user to show different channel settings. If ALL is selected, then all four-channel settings are displayed at once.</p> <p>Note: The following settings, denoted by Ch X, are the same for all four channels.</p>
Ch X Power Limit	<p>Options: On, Off</p> <p>This button controls the individual channel. This allows the user to turn off an individual channel, even if the Output Power is on.</p>
Ch X Power Limit Slider and Numeric Up Down	<p>Range: 0-100.0</p> <p>This sets the power limit of an individual channel. This scaler is applied to the main power setting to calculate the actual power of a channel.</p> <p>Note: This setting can be lowered to limit the total output power. This is useful in situations where user control of the front knob is required, but a maximum light output is needed at less than 100%.</p> <p>Note: The default setting on single channel units is 75%, to limit temperature of the LED while at minimum fan speed. This can be increased by changing to Industrial or Heavy Industrial mode, as outlined in the Unit Operating Modes section on page 45.</p>
Ch X Shutdown Pin	<p>Options: Active Low, Active High</p> <p>This setting controls when the unit will be in shutdown mode based off the input from the multiport. The default is Active Low, so if the Multiport Shutdown Pin is pulled to GND (either driven or with a contact closure), the</p>

LED channel will not turn on. In a single channel device, a low signal on Channel 1 will turn off the light completely.

Continuous Strobe

In Continuous Strobe mode, the unit will generate a trigger source from its internal clock. This allows the user to set up a continuous strobe train to stop motion in machine vision applications.

Note: Strobe intensity is set by the analog dimming level under the “LED Channel Settings” section.

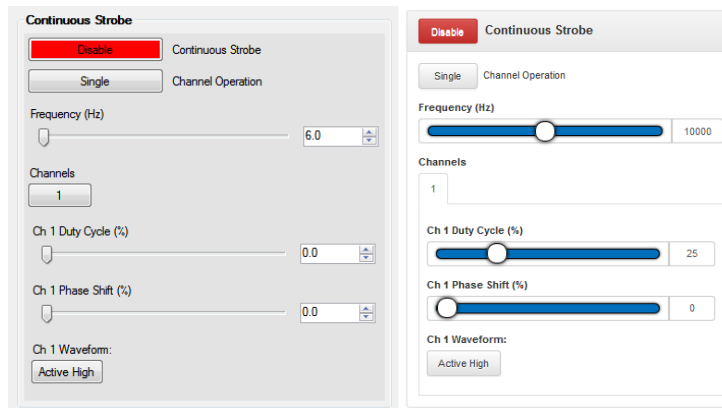
Figure 10: Continuous Strobe (Left: Dashboard, Right: Web Interface)

Control Name	Description
Continuous Strobe	Options: Enable, Disable When this option, along with the Output Power, is enabled, the LED will strobe at the settings listed below.
Channel Operation	Options: Single, Quad When in Single mode, all four channels strobe with the settings of Channel 1. Quad mode shows settings for each channel individually.
Frequency Slider and Numeric Up Down	Range: 6 – 20,000 Hz Frequency of the strobe trigger.
Channels	Options: 1, 2, 3, 4, All When in Quad mode, all options are displayed to allow the user to show different channel settings. If ALL is selected, then all four channel settings are displayed at once. Note: The following settings, denoted by Ch X , are the same for all four channels.
Ch X Duty Cycle (%)	Range: 0-100.0 Percent of the waveform that is active.
Ch X Phase Shift (%)	Range: 0-100.0 Percent of the waveform the strobe will wait before becoming active.
Ch X Waveform	Options: Active High, Active Low

Sets the state of the LED during the Active part of the waveform. High = LED On, Low = LED Off.

Below are two samples of Continuous Strobe trains. The LED output is detected on a reverse biased photodiode. In Figure 11 below, because the Waveform is set to “Active High”, the LED is ON for only 25% of the waveform. In Figure 12 below, the Waveform is set to “Active Low”, the LED is OFF for only 25% of the waveform.

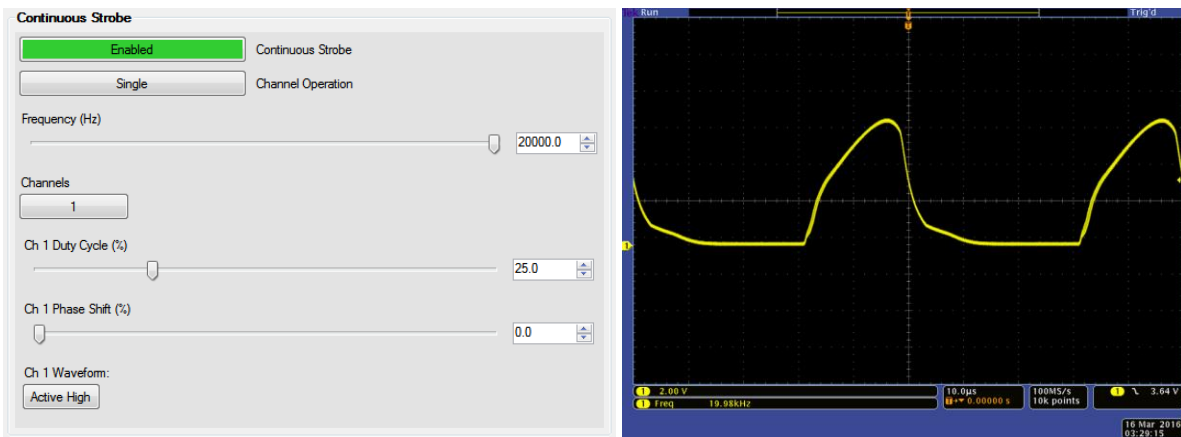
Figure 11: Active High Continuous Strobe Train (Left: Strobe Settings, Right: Light Output measured with Photodiode)

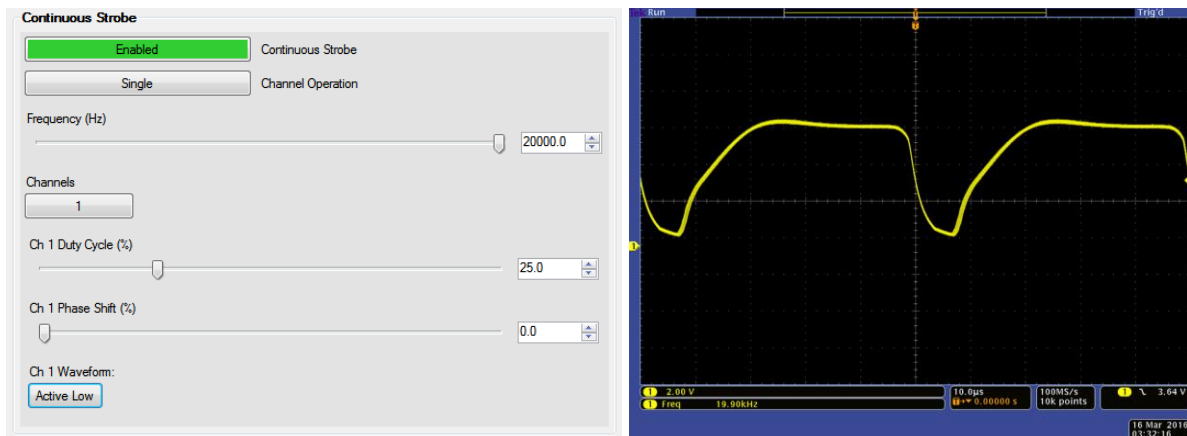
Figure 12: Active Low Continuous Strobe Train (Left: Strobe Settings, Right: Light Output measured with Photodiode)

Triggered Strobe

In Triggered Strobe mode, the unit will wait for a user generated trigger signal on the digital pins of the multiport. This allows the user to set up a strobe train, synced with an external trigger source, to stop motion in machine vision applications.

Note: Strobe intensity is set by the analog dimming level set under “LED Channel Settings” section.

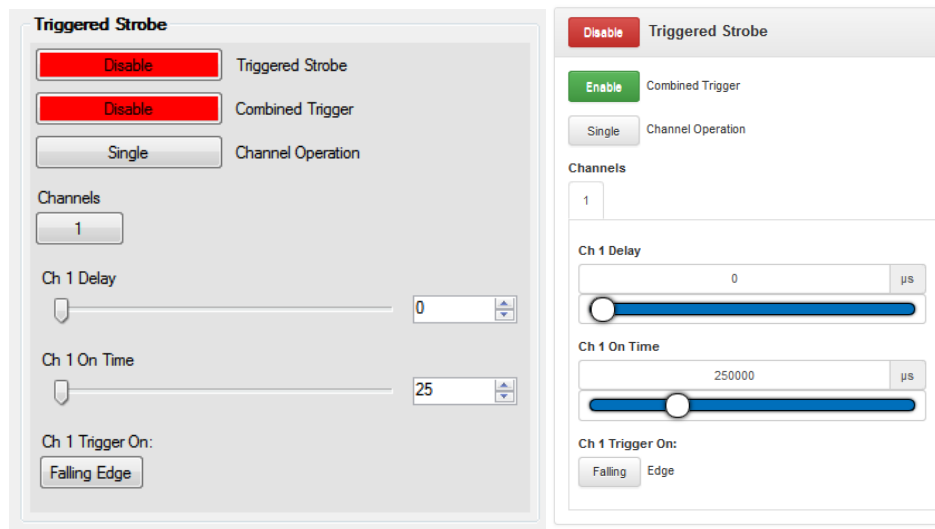
Figure 13: Triggered Strobe (Left: Dashboard, Right: Web Interface)

Below is a sample strobe train. The input is a 20 KHz, 0V Low, 3.3V High signal. The output is detected on a reverse biased photodiode. The spread on the output pulse train is caused by the input signal falling at different times during the 5μs tick of the internal clock used to control the LED.

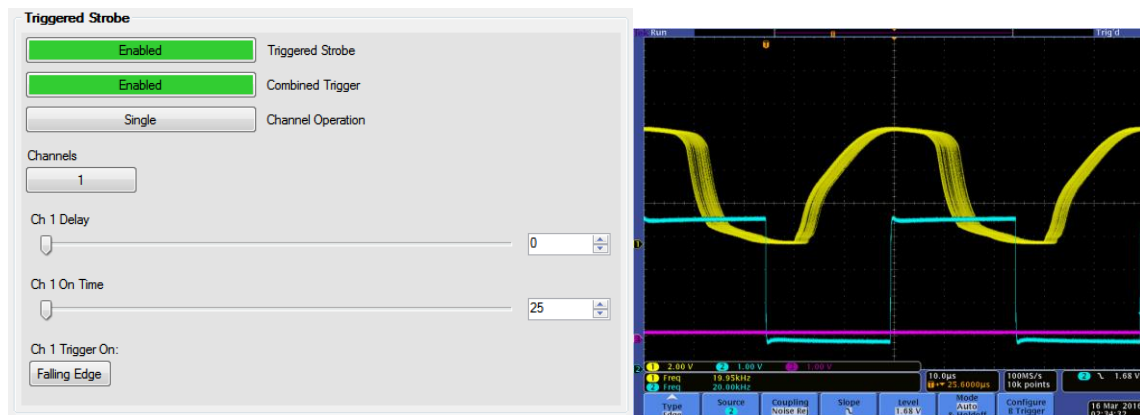
Figure 14: Triggered Strobe Train (Left: Strobe Settings, Right: Oscilloscope trace, Bottom 20 KHz input, Top Light Output measured with Photodiode)

Control Name	Description
Triggered Strobe	Options: Enable, Disable When this option, along with the Output Power, is enabled, the LED will strobe at the settings listed below.
Combined Trigger	Options: Enable, Disable When this option is enabled, the system will trigger a strobe on all channels when there is a signal on any of the four multi-port digital inputs.
Channel Operation	Options: Single, Quad All four channels strobe with the settings of Channel 1 when in Single mode. Quad mode shows settings for each channel individually.
Channels	Options: 1, 2, 3, 4, All When in Quad mode, all options are displayed to allow the user to show different channel settings. If ALL is selected, then all four channel settings are displayed at once. Note: The following settings, denoted by Ch X , are the same for all four channels.
Ch X Delay Slider and Numeric Up Down	Range: 0 – 1000000µs Number of microseconds the channel will delay, from when the trigger signal is received, before turning on the LED. Note: This setting is in increments of 5µs. You can set a value in increments of 1µs for Legacy Command response, but the system will round down to the nearest 5µs increment. Note: There is a 4 – 9µs delay in the response time of the system. This setting adds to that time.
Ch X On Time Slider and Numeric Up Down	Range: 25 – 1000000µs Number of microseconds the channel will remain on for the strobe pulse. Note: This setting is in increments of 5µs. You can set a value in increments of 1µs for legacy command response, but the system will round down to the nearest 5µs increment.
Ch X Trigger On	Options: Falling Edge, Rising Edge Channel will trigger on the falling or rising edge of the input signal, on the digital pin of the multiport, depending on this setting.

Equalizer

The CV-LS Equalizer is a closed loop controller using a small light tap off the led to control the source current and maintain a constant power output. This value is displayed as the “Light Feed Back” (LFB) in the Status section, and “Current Light Output” (CLO) in the Equalizer section. The LFB value is calculated from a much shorter averaging period than the CLO value. The LFB can be used to see a live view of the light output, while the CLO value is used by the internal control loop of the Equalizer.

The basic operating modes of the light source use constant current supply for the LED. This maintains a constant output over time as long as the LED temperature does not change. Age also plays a role, but in LEDs, the timespan is quite long. The LED in the CV-LS is rated for 50,000 hours at an LED temperature of 70°C. This means that at 50,000 hours the LED will still produce 70% of its intensity.

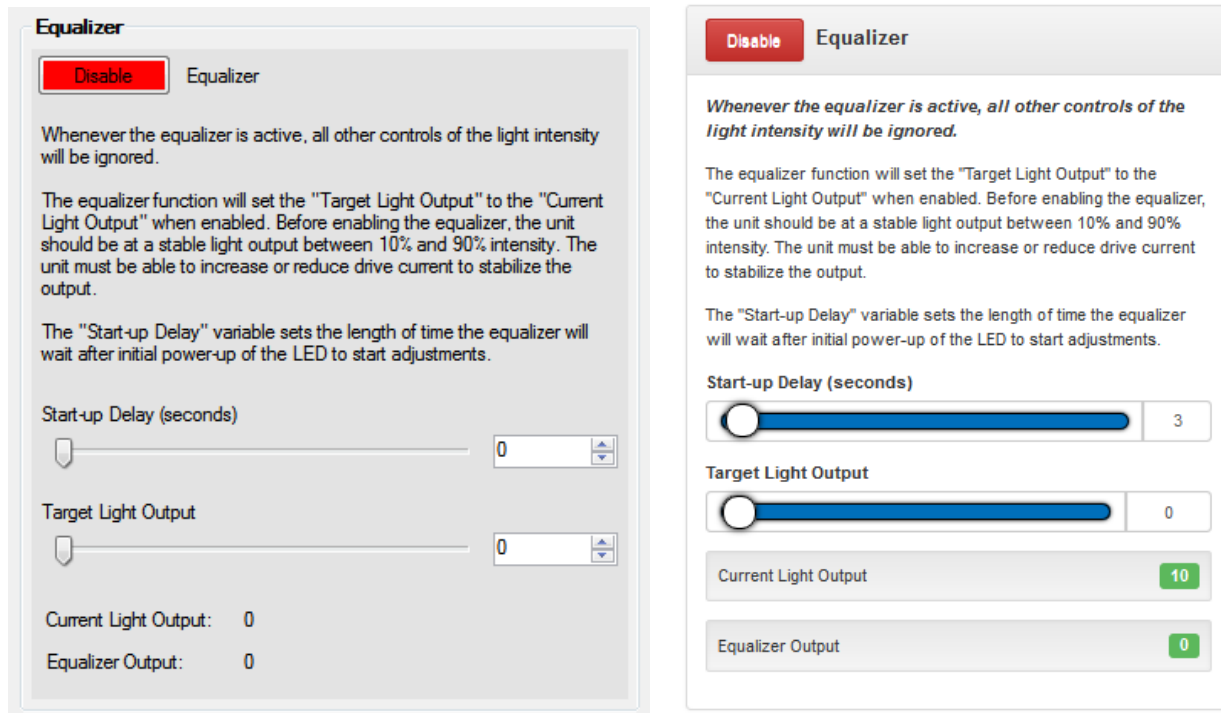
Figure 15: Equalizer (Left: Dashboard, Right: Web Interface)

Equalizer Key Points

Using the equalizer has limitations, which should be understood before using it.

First, the equalizer relies on the CLO, which is averaged over many milliseconds to maintain a steady control loop. Because of this, the equalizer may not perform well if used with strobe modes.

Second, the equalizer needs enough room in the driver to increase or decrease the current with time to maintain the light output. This means if the light is already at 100%, the equalizer cannot increase the current if light output decreases.

Third, because the light output will change significantly when the unit is first turned on, the user should give sufficient time for the light output to stabilize at a satisfactory level before engaging the equalizer. When the equalizer is enabled for the first time, it will set the current light output as its target. This can be manually changed later with the “Target Light Output” controls if needed.

Once the settings are stored, upon future power ups, the unit will return to the set “Target Light Output” value.

The final control in this section is to a time delay. This is a delay, in seconds, after the unit “Output Power” is enabled, before the equalizer will start controlling the output current.

Fan Control

The CV-LS has a variable speed fan to optimize fan noise versus cooling needs. The higher fan speeds allow for operating the LED at max intensity under warmer ambient temperature, but is much louder. The minimum fan speed still allows 70% light output at room temperature and is much quieter.

The user can test the noise of the fan by changing the “Fan Control” from “Auto” to “Manual”. The manual mode will allow for the fan to be on at the speed shown, even if the LED is off. The fan will stay in “Manual” mode for 15 seconds after each speed change, and will then change back to “Auto” mode.

The status windows can be used to monitor the LED temperature to find the optimal tradeoff of fan speed and LED temperature for the intended application.

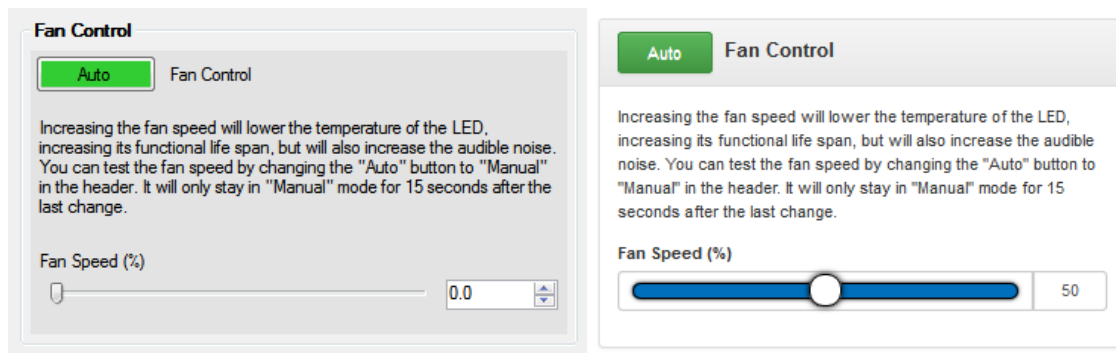
Figure 16: Fan Control (Left: Dashboard, Right: Web Interface)

Upgrade (ADMIN Level)

The Upgrade sections allow the user to update to the latest firmware without sending the unit in under an RMA. Upgrading the firmware will not change any of the unit settings, or disrupt the settings stored at the factory. If upgrading a unit causes any issues, please report the problem to technical support for evaluation.

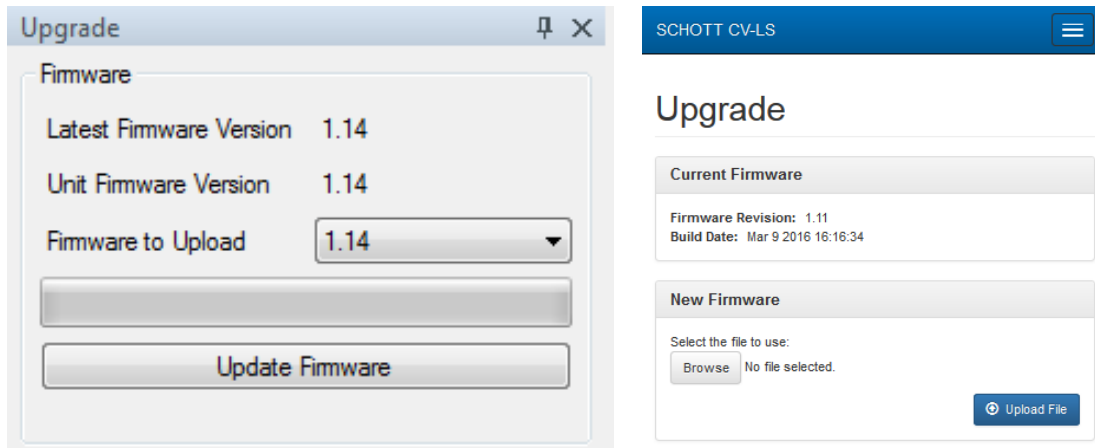
Figure 17: Upgrade (Left: Dashboard, Right: Web Interface)

Control Name	Description
Update Firmware	The “Latest Firmware Version” is the latest firmware built into the Dashboard. The user can select which firmware version to upload by changing the value in the drop list “Firmware to Upload”. The user can manually upgrade the connected unit to the selected firmware by clicking the “Update Firmware” button.
Upload File	In the Web Interface, clicking this button will open a file dialog box, allowing the user to select a firmware .bin file to upload to the unit. This can be used to downgrade a firmware if required.

Settings (ADMIN Level)

The Settings section of the controls are static pages on all interfaces. Any changes that are made on this window must be saved to memory before they can be made live before they take effect. This is because the settings on this page require restarting core parts of the processor, and cannot be done without disrupting the connection to the user.

The following sections will describe in detail the controls in each part of the Settings windows.

Note: To save changes, click “Write All” in the Dashboard, or “Save” in the Web Interface.

Note: The controls are nested inside tabs in the Web Interface, but are tiled down the page in the Dashboard Window.

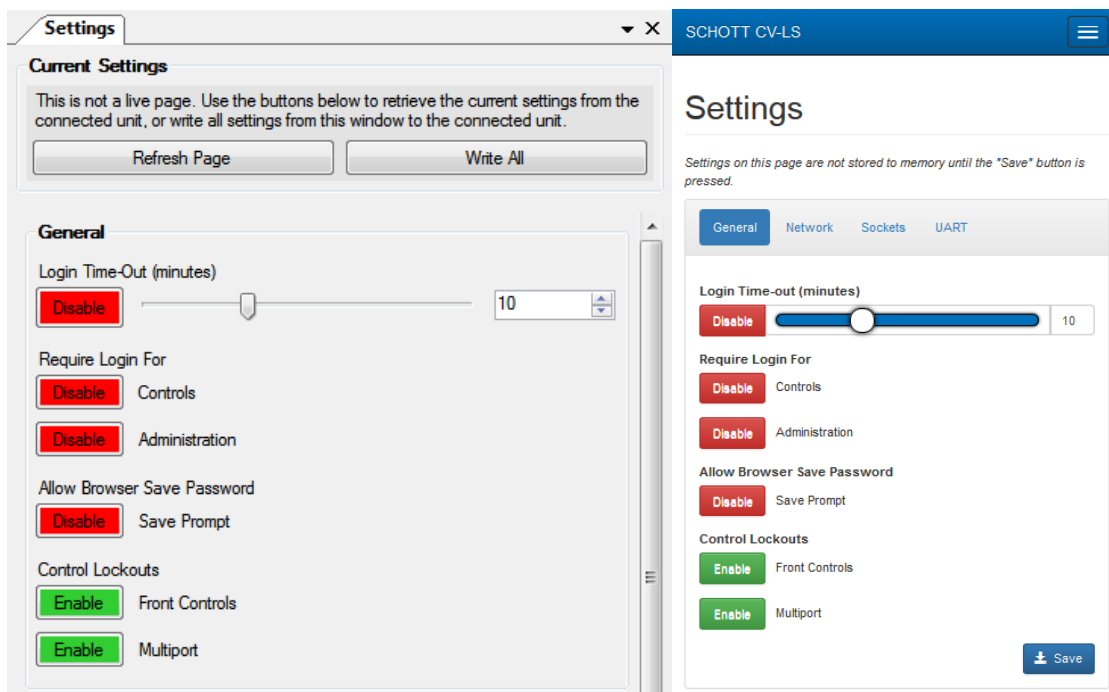
Figure 18: Settings (Left: Dashboard, Right: Web Interface)

General

The General section controls how users log into the device, and which interfaces are active.

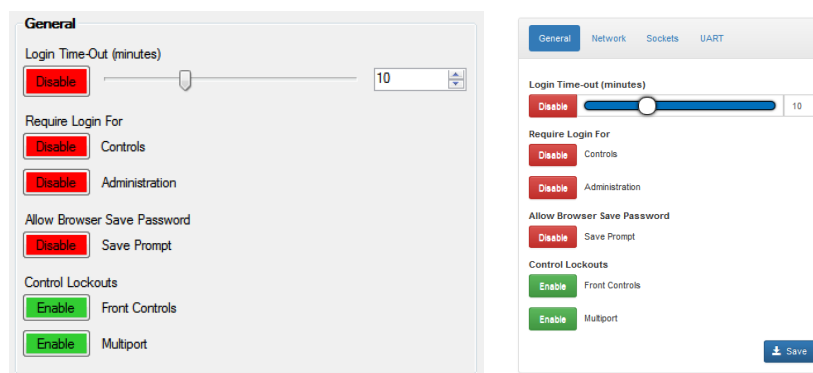
Figure 19: General (Left: Dashboard, Right: Web Interface)

Control Name	Description
Login Time-Out Enable	Options: Enable, Disable When enabled, the user will only stay logged in for a pre-set time before the unit will log them out.
Login Time-Out Slider and Numeric Up Down	Range: 1 – 30 minutes If logins are enabled, this time is the delay before the user is automatically logged out. Anytime a new page is viewed or a setting is changed, this timeout is renewed to the full time. Note: The user can also manually log out using the controls available.
Require Login for Controls	Options: Enable, Disable When enabled, the system will require a login of USER level or higher to use USER pages. The Controls page is currently the only page requiring a USER level login.
Require Login for Admin	Options: Enable, Disable When enabled, the system will require a login of ADMIN level to use ADMIN pages. In the Web Interface, any pages under the Administration menu require an ADMIN level login. These include: Settings, Configurations, Users, Upgrade, and Diagnostics.
Allow Browser Save Password	Options: Enable, Disable When enabled and a user logs into the Web Interface, the browser will prompt to save the password.
Lockout Front Controls	Options: Enable, Disable When enabled, the unit will ignore the front button and knob, except when using the front button to performing a factory reset.
Lockout Multiport	Options: Enable, Disable When enabled, the unit will ignore analog signals to the multiport. Digital signals are still processed.

Network

The Network section allows the user to specify how the unit will connect to a network. By default the units have the following settings:

- Hostname: SCHOTT CV-LS
- DHCP: Enabled
- IP Address: 192.168.0.2
- Subnet Mask: 255.255.255.0
- Default Gateway: 192.168.0.1
- Primary DNS: 192.168.0.1
- Secondary DNS: 192.168.0.1

This will allow the unit to work without changing these settings on most networks. If these settings are not compatible with your network, it is easiest to change the network settings using the Dashboard connected by USB.

A useful feature of the network pages are the buttons to the right of each field. These buttons display the current network settings, and copy their contents into the field on the left when clicked. This allows you to store the current DHCP settings, and then disable the DHCP to have a static configuration with the values assigned by the DHCP server on your network.

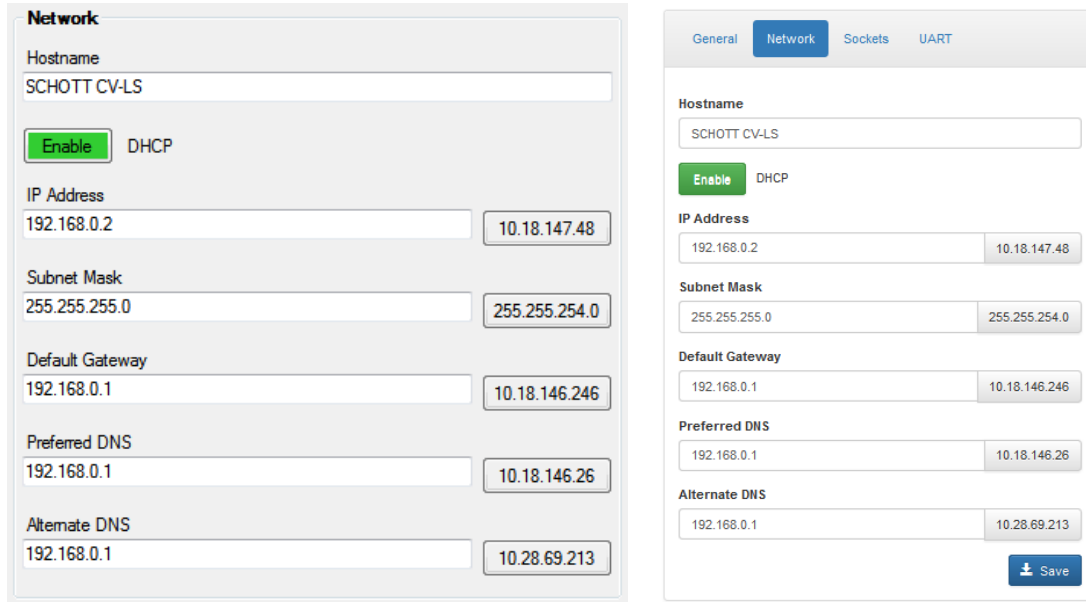
Figure 20: Network (Left: Dashboard, Right: Web Interface)

Control Name	Description
Hostname	The name that will display in the finder application. Used to easily identify multiple units on the same network.
DHCP	<p>Options: Enable, Disable</p> <p>When enabled, the unit will first request network information from the local DHCP server. If this fails, it will fall back to the static network information stored in memory.</p> <p>Note: The unit will be available more quickly on the network if DCHP is disabled when no DHCP server is present. Otherwise, a timeout must occur before falling back to the static IP in memory.</p>
Text Box Values	The static network values of the system.
Button Text Values	<p>The current DHCP values assigned to the unit.</p> <p>Note: These will be 0.0.0.0 if DHCP is disabled.</p>

Sockets

The sockets are used to communicate with the unit over the network programmatically. The Legacy Socket is a standard text based telnet socket, and uses the same command structure as the UART and USB interfaces.

The Binary Ethernet Socket is a new interface based on a binary protocol with packets and error checksums. This interface is much faster and easier to program, and is best used for continuous interaction with the unit.

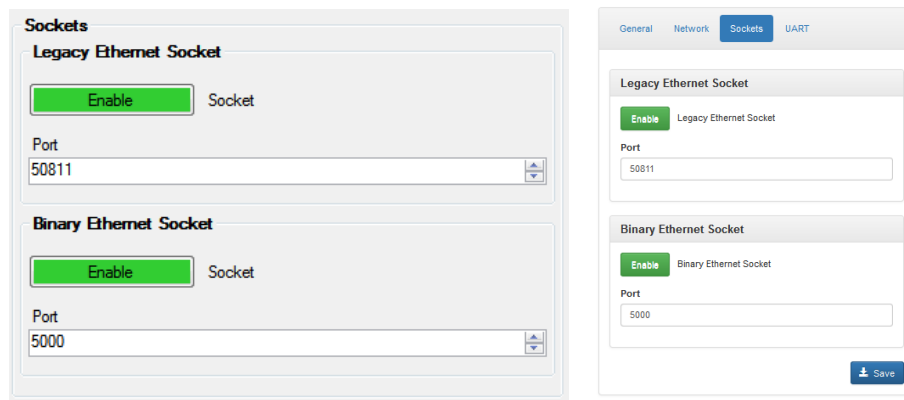
Figure 21: Sockets (Left: Dashboard, Right: Web Interface)

Control Name	Description
Socket	Options: Enable, Disable Port is only active when this setting is Enabled. Unit will refuse connections when Disabled.
Port	Range: 1 - 65,536 The port of one socket must be different than the port of the other socket.

UART

The UART settings allow the user to change the communications protocol of the Legacy UART pins in the Multiport. The default settings are the most common UART port and should only be changed if the user has purposely changed the settings in their command terminal.

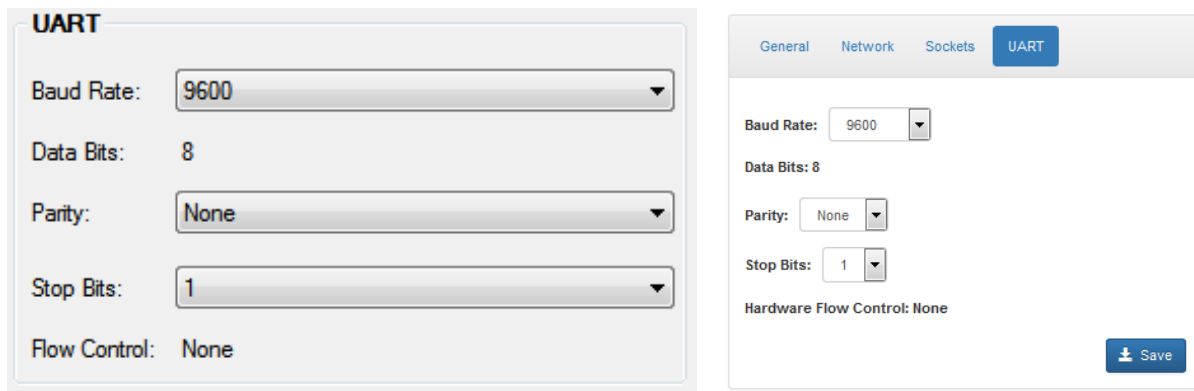
Figure 22: UART (Left: Dashboard, Right: Web Interface)

Control Name	Description
Baud Rate	This is the communications speed. The higher the number, the faster the unit will be able to receive and transmit text. However, not all computers / UART adaptors can run at the highest speed. If you cannot connect at a higher speed, work up to the highest speed your computer can use.
Parity / Stop Bits	These settings rarely need to be changed. They affect the protocol used to transmit characters.

Users (ADMIN Level)

The CV-LS unit comes with two pre-set users, shown in Figure 23 below. Each unit can have up to 10 users.

- 1) admin
 - a) User Name: admin
 - b) User Password: admin
 - c) Account Level: ADMIN
- 2) user
 - a) User Name: user
 - b) User Password: user
 - c) Account Level: USER

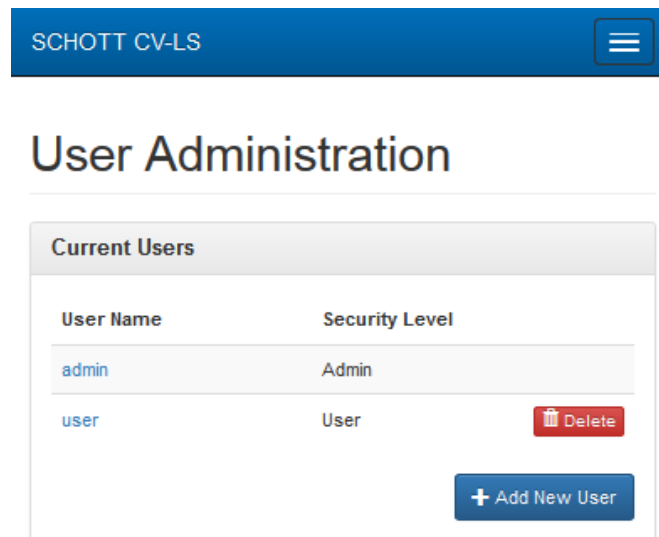
Figure 23: Users

These accounts can be modified as needed for the intended application by clicking on the “User Name” column. There must always be an ADMIN account in the first spot, and this cannot be removed. Any other account can be deleted with the “Delete” button to the right of the user.

New users can be added with the “Add New User” button at the bottom of the screen.

Add / Edit User

When adding or editing a user, different fields are required. If any field is incorrect, a message will be displayed with the error.

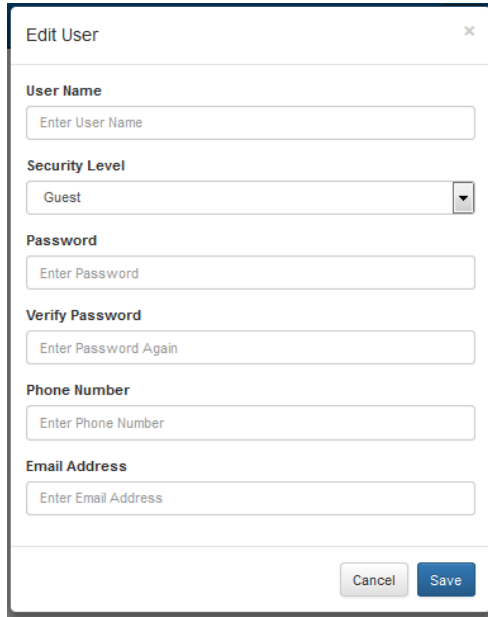
Figure 24: Add / Edit User

Control Name	Description
User Name	Required For: Add, Edit This used to log in.
Security Level	Required For: Add, Edit This limits what the user can view when logged in.
Password / Verify Password	Required For: Add When editing a user, leaving this field blank will keep the old password.
Phone Number	Required For: None This field is for a planned firmware upgrade, where software errors / exceptions will cause a notification to be texted to the ADMIN level phone numbers stored in the unit.
Email Address	Required For: None This field is for a planned firmware upgrade, where software errors / exceptions will cause a notification to be emailed to the ADMIN level emails stored in the unit.

Delete User

When clicking the “Delete” button to the right of the user, the confirmation dialog shown in Figure 25 below will be displayed. Once the “Delete” button on this dialog is clicked, the user will be deleted and non-retrievable.

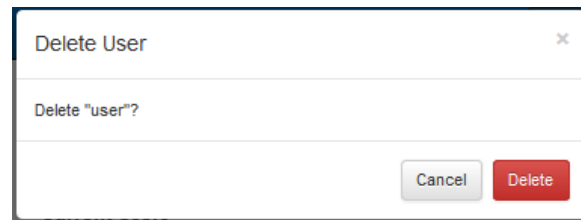
Figure 25: Delete User

Diagnostics (ADMIN Level)

The Diagnostics window is used when something unexplainable is happening to the unit. Information on this page will help technical support diagnose any issues you are having with a unit.

The Logs section has diagnostic information showing where the code was when the exception occurred. This information will allow tech support to track down any bugs in the firmware. The log can be cleared or exported to a text file with the buttons below the table.

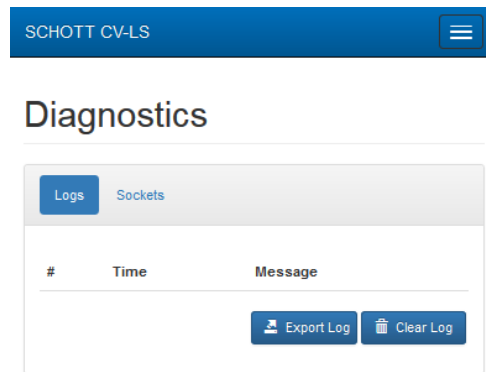
Figure 26: Diagnostics

Sockets

The Sockets tab will show if a computer is currently connected to one of the sockets, and allow an ADMIN level account to kick that computer off the socket. This is useful if someone left the Dashboard connected to a unit remotely and a different user needs to access it. That user can then use the web interface to kick them off the socket, and connect to the socket themselves.

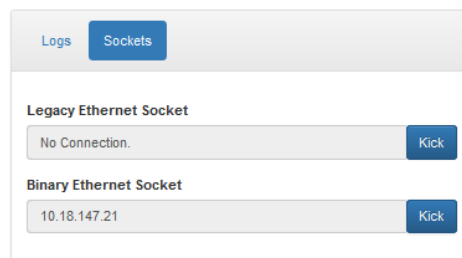



Figure 27: Sockets

Legacy Socket

Legacy Socket

Socket Connection

The Legacy Socket can be connected to with any telnet type application. TeraTerm is one such application. Once the IP address of the unit is known, follow the steps below to connect to the Legacy Socket in TeraTerm.

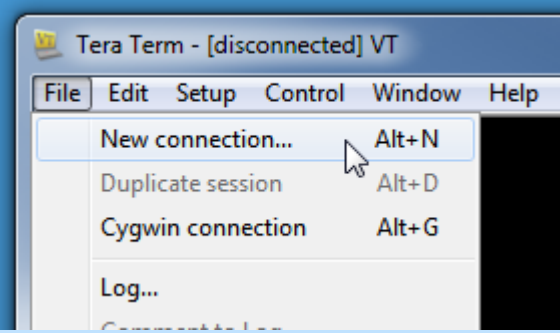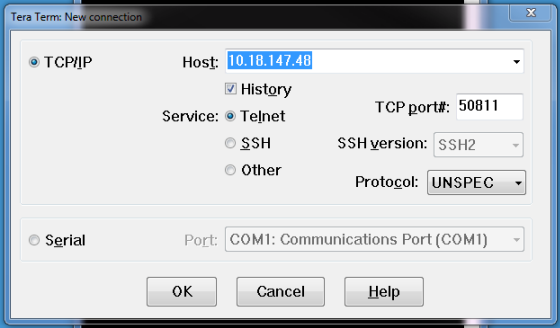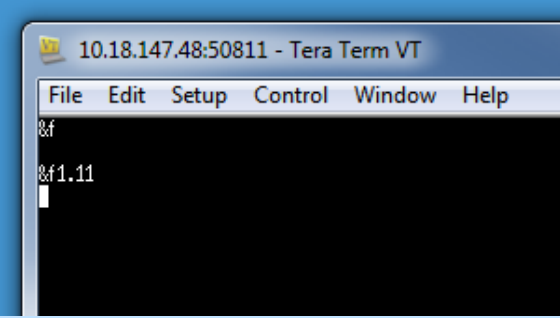Steps	Description
	With TeraTerm open, start a “New Connection”.
	<ol style="list-style-type: none">1. Enter the IP address of the unit into the “Host” field.2. Enter the port for the Legacy Socket into the “TCP port#” field (Default is 50811).3. Click “OK”.
	<p>The title of the window should change to the IP address and port number entered in the previous step. This means TeraTerm is connected to the socket.</p> <p>Legacy commands can be typed to control the unit. The “&f\r” command is shown here to retrieve the firmware version number (V1.11 in this case).</p>

Command Structure

All commands start with the ampersand character “&” and end with a carriage return “\r”. There is a string representing the command immediately after the “&”, followed by the desired value.

The unit will discard all characters until it receives an “&” start character. Once the carriage return “\r” is received, the unit will process the command. If a valid command is not found, the unit will return “&n” followed by the command string with an indicator of which character was not recognized.

Negative Acknowledgements

Negative acknowledgements are used to indicate that a received command was invalid. The following negative acknowledgements are supported:

Invalid Command:	Return string format: &nXpY &n = Negative acknowledgement X = Received characters parsed correctly p = Invalid character location start Y = Invalid character received
Invalid Parameter:	Return string format: &nXpYYY &n = Negative acknowledgement X = Received characters parsed correctly p = Invalid parameter location start YYY = Invalid parameter received

Commands

Please download the programming guide found under the Downloads Section on page 94 for a full list of available commands and how to use them.

Binary Socket

Socket Connection

The Binary Socket can be connected to using custom applications that allow for fast and robust communications. The Dashboard Application uses the binary socket to enable live views and graphing at high speed. The easiest way to start working with the Binary Socket is to look at the source code of the Dashboard Application. This section of the manual describes some of the core functions used for this interface, but is not meant as a sole source of information from which to create an interface.

Note: Example code is copied from the Dashboard source code, and is written in C#.net.

Note: Please reference the source code of the Dashboard Application and the Programming Guide for supplemental information to create your own implementation of the Binary Socket.

Fletcher16 Checksums

The Fletcher16 Checksum is used to determine if a packet is corrupted or not. There are many ways to calculate the Fletcher16 checksum, but this is the implementation used in the Dashboard Application:

```
// computes the fletcher16 checksum for the data[]
public static byte[] fletcher16(byte[] data, int bytes)
{
    UInt16 sum1 = 0xff, sum2 = 0xff;
    int tlen;
    int i = 0;

    while (bytes > 0)
    {
        tlen = bytes > 20 ? 20 : bytes;
        bytes -= tlen;
        do
        {
            sum2 += sum1 += data[i++];
        } while (--tlen > 0);
        /* First reduction step to reduce sums to 8 bits */
        sum1 = (UInt16)((sum1 & 0xff) + (sum1 >> 8));
        sum2 = (UInt16)((sum2 & 0xff) + (sum2 >> 8));
    }

    /* Second reduction step to reduce sums to 8 bits */
    sum1 = (UInt16)((sum1 & 0xff) + (sum1 >> 8));
    sum2 = (UInt16)((sum2 & 0xff) + (sum2 >> 8));

    return new byte[2] { (byte)sum2, (byte)sum1 };
}
```

Encoding a Packet

All binary commands have the following key Byte Values:

Byte Name	Value
Start	0x12
Escape	0x7D
End	0x13

A command has the following structure:

- 1) 2 Byte Command
 - a) Bits 1-12: Command Code
 - b) Bit 13: Read when 0, Write when 1
 - c) Bits 14-15: Command Set
- 2) 2 Byte Data Length (number of bytes in Data Payload)
- 3) Data Payload
- 4) 2 Byte Fletcher16 Checksum

Once the command is constructed, the data must be escaped. Any byte in the entire packet that matches the **Start**, **Escape**, or **End** bytes must have the **Escape** byte added just before it. The byte must then have a bitwise XOR with 0x80 applied. This ensures that no byte in the entire payload will match the **Start** or **End** bytes.

Once the command is escaped, a **Start** byte is prepended to the packet and a **Stop** byte is appended to the packet.

This process creates a single complete packet that can be transmitted to the Binary Socket. If a packet does not follow this form, it will be discarded by the unit.

Note: An example implementation of this method can be found in the “classBinarySocket.c” file in the Dashboard source code.

Decoding a Packet

Decoding a stream of data from the unit can be done by reversing the process of “Encoding a Packet”. The Dashboard Application implements an asynchronous approach to decoding packets. This is the preferred method to make use of the streaming status functions available to the Binary Socket.

Commands

Please download the programming guide found under the Downloads Section on page 94 for a full list of available commands and how to use them.

Product Specification

Environment

Operational Temperature

CV-LS: 0°C to 45°C.

Power Supply: -30°C to 50°C.

There are temperature sensors on the LED module and PCB. Readings are accessible through software interfaces. Software controls will not allow the LED to be driven over its maximum operating temperature and will automatically dim the LED to prevent damage.

Storage Temperature

CV-LS: -25°C to 85°C

Power Supply: -40°C to 85°C

Humidity

CV-LS: 5% to 95% Non-Condensing

Power Supply: 20% to 90% Non-Condensing

Electrical

Input

CV-LS: 18-28VDC / 5A Max

Power Supply: 85 – 264 VAC, 47 – 63 Hz, 120W

CAUTION

The CV-LS must have an Earth Ground connection for EMC Compliance.

Users must adhere to one of the following:

- Use the included AC/DC power supply. Third prong must be connected to Earth Ground.
- One of the GND pins in the Multiport Connector must be connected to Earth Ground.

Redundant Power Inputs

The CV-LS can be powered through either the Multiport terminal block or the DC Power Input. Both of these inputs are isolated from the other through a load sharing circuit built into the CV-LS. This means the user can supply power through both interfaces without damaging either supply, allowing for redundant supply inputs.

Product Specification

Mechanical

Weight (Typical)

CV-LS: 1.8 lbs

Power Supply: 1.7 lbs

Packing Materials: 1 lbs

Total Shipping Weight: 4.5 lbs

Physical Dimensions (Inches)

The following diagram shows the physical dimensions of the CV-LS.

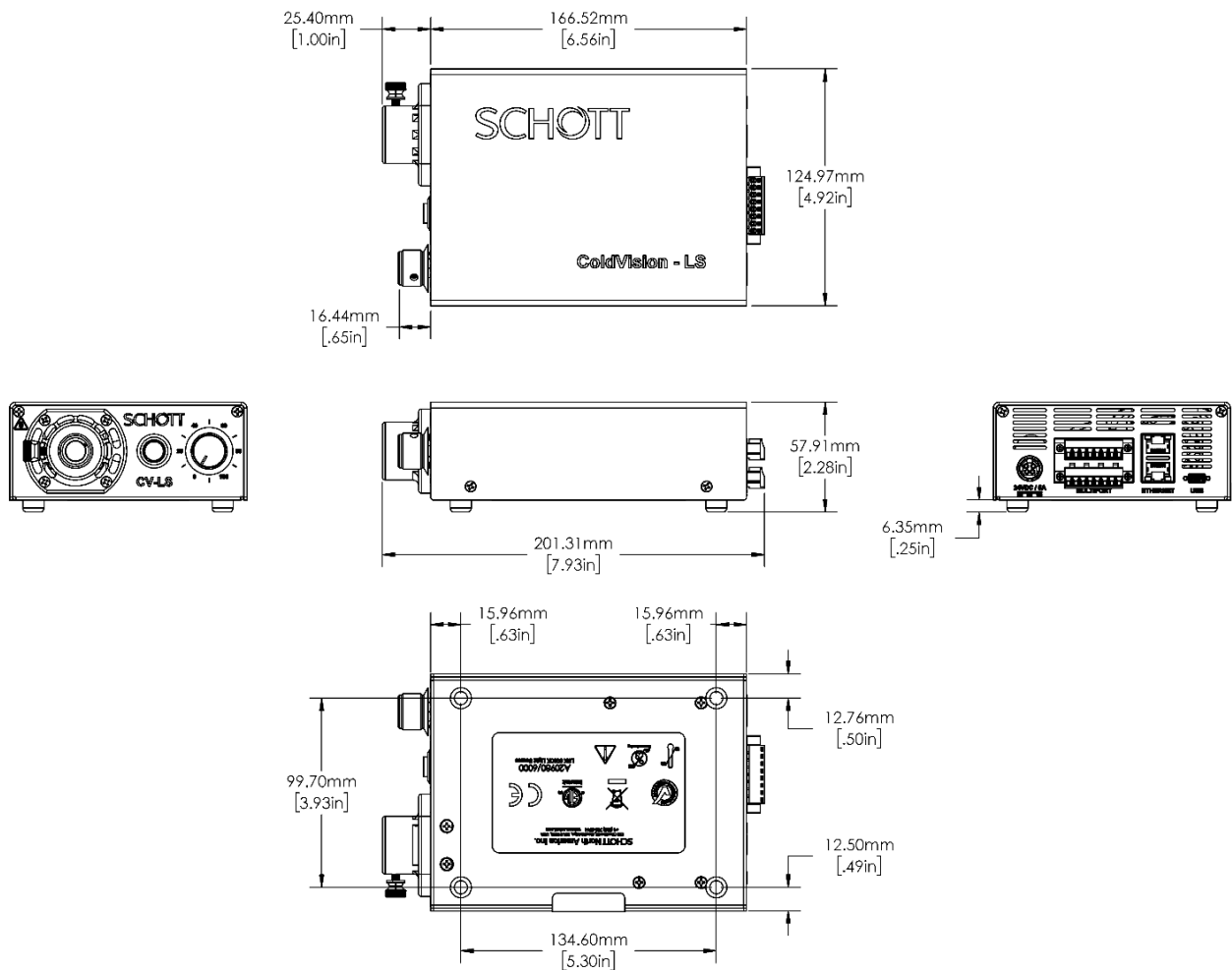
Figure 28: CV-LS Dimensions

Product Specification

Front Panel Description

The front of the unit is where the manual controls for use by an operator exist. This is also the exit point for the light, as well as the ventilation direction for air exiting the LED heat sink. These vents should never be impeded, as it may result in reduced light output.

CAUTION

The nose cone and surrounding area may become hot at high light output. Under certain user settings, these surfaces may become hot enough to burn the operator. Always handle these surfaces accordingly.

Figure 29: CV-LS Front Panel Description

Label	Description
Vent Outputs	Heated air from the LED will vent out of these slots. They must remain unobstructed to maintain full light output.
Fiber Optic Thumb Screw	This thumbscrew is used to retain fiber optic cables inserted into the nose cone.
Output Control Knob	This knob allows for overall adjustment of the LED light output.
Power Switch / Indicator	This switch can be used to enable/disable the LED on the unit. The LED indicator ring will turn on when the LED is enabled, or flash if there is an error with the unit.

Product Specification

Rear Panel Description

The back of the unit is where power is supplied, all remote control interfaces are connected, and the air intake vents for LED cooling are located. All ports on the back of this unit are locking connectors for industrial use.

CAUTION

The vent slots on the rear panel supply air to cool the LED. Ensure adequate airflow to vents at all times.

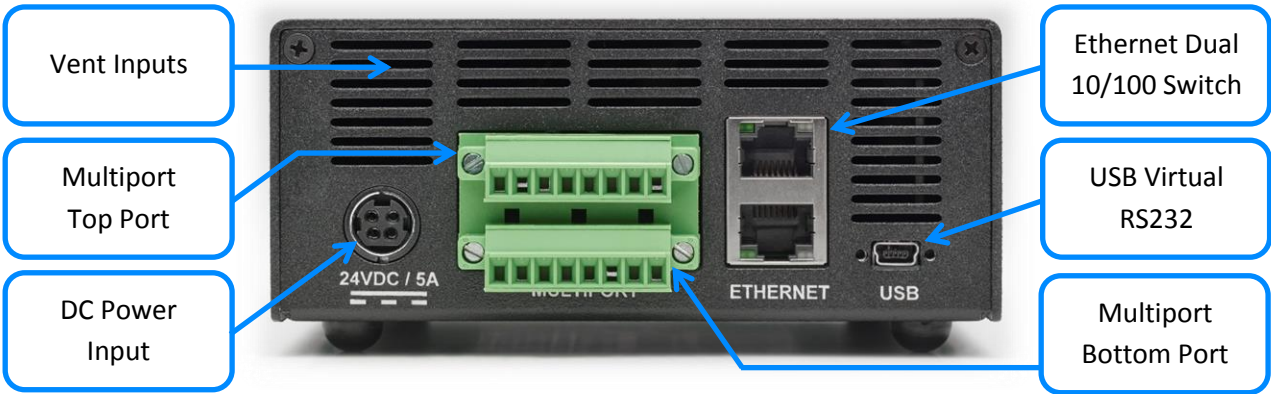
Figure 30: CV-LS Rear Panel Description

Label	Description
Vent Inputs	Air is pulled into these vents to cool the LED.
DC Power Input	The included AC/DC power supply will plug into this locking connector.
Ethernet Dual 10/100 Switch	The dual Ethernet port is a built in 10/100 Ethernet Switch. Either port can be used to connect the unit to a network. The second port can be used to chain units off the same network port.
USB Virtual RS232	This port will be detected as a RS232 port and can be used in any software that would normally connect to RS232. (See the Downloads section on page 94)
Multiport Top Port	This connector contains CH 2 to CH 4 analog and digital inputs, as well as the native RS232 TX and RX pins.
Multiport Bottom Port	This connector contains CH 1 analog and digital inputs, fault output, DC power input, 5V output, and GND connections.

Multiport Pin Details

⚠ CAUTION

All wiring to the Multiport terminal block signal lines must be rated to 30V/200mA. The +24V and GND wiring must be rated for at least 50V/5A. The temperature rating of all wiring should be greater than 70°C. The Multiport terminal block is rated for 16-28 AWG solid or stranded wire.

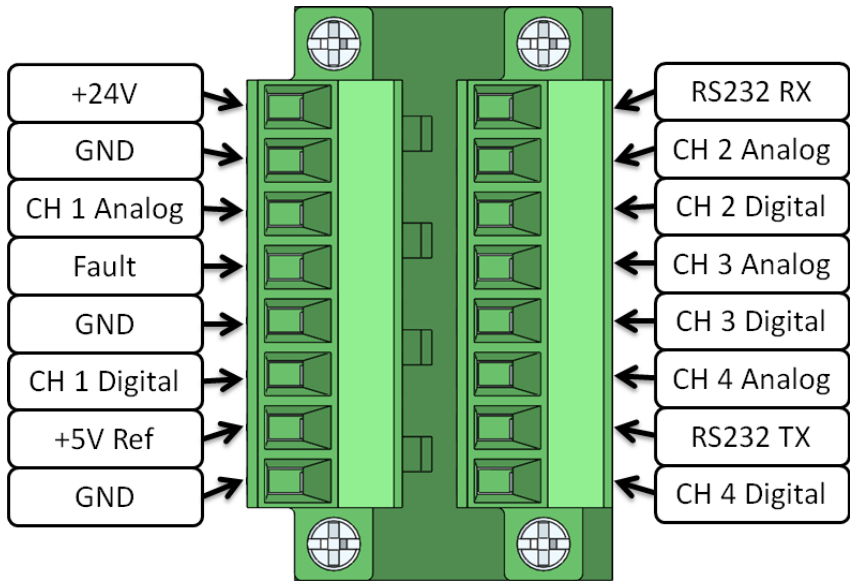
Figure 31: CV-LS Multiport Description

Name	Voltage Range	Notes
+24V	18V to 28V	Power connection must be a low impedance source. Current ramp rate can be as much as 4A in 25µs.
GND	Earth Ground	When powering the unit through the multiport, at least one GND pin should be connected to Earth Ground through a low impedance path.
CH X Analog	0V to 5V	Controls the intensity of the LED channels.
CH X Digital	0V to 24V (3.3V TTL, 24V Max)	Pin is pulled up internally; a contact closure can be used to create a low signal. <0.5V = Low
Fault	Open Collector (24V Max)	Active Low: Signal will be pulled low (GND) when a fault exists.
+5V Ref	5V Output Reference	Voltage rail can be used to power potentiometers for analog inputs.
RS232 RX	-30V to +30V	Data input to the device.
RS232 TX	-5V to +5V	Data output from the device.

LED Information

 **CAUTION**

Risk Group 2

CAUTION: Possibly hazardous optical radiation emitted from this product.
Do not stare at operating lamp.
May be harmful to the eyes.

Light output is measured using a SCHOTT fiber optic light guide, Ø 13 mm active, length 1 m (A08051.40 bundle).

LED Light Output will vary slightly with LED temperature. Please allow time for the temperature to equalize to achieve stable light output. This can take several minutes at maximum fan speed settings.

LED output is through the front nosepiece of the unit. Please follow all safety and operational guidelines to ensure safe operation of the unit.

A20980/6000K

Typical Output: 1300 lumens.

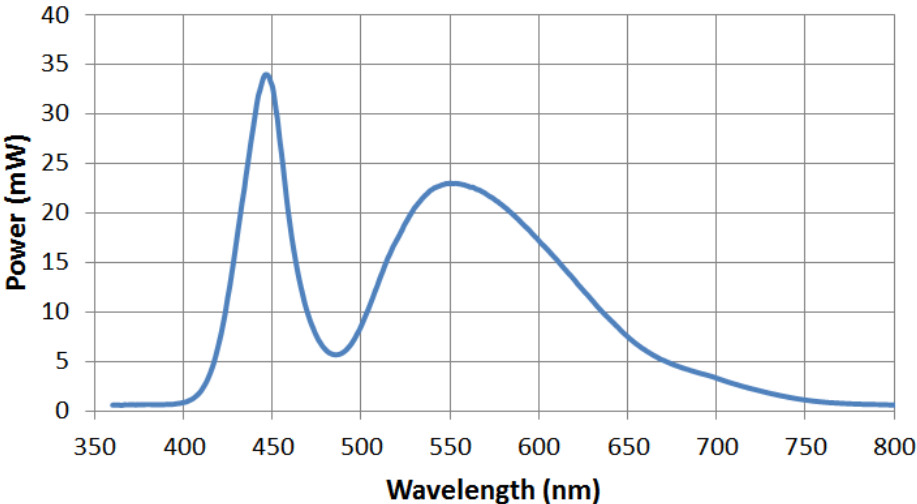
Figure 32: Typical relative spectral power vs wavelength (6000K model)

A20980/3000K

Typical Output: 900 lumens.

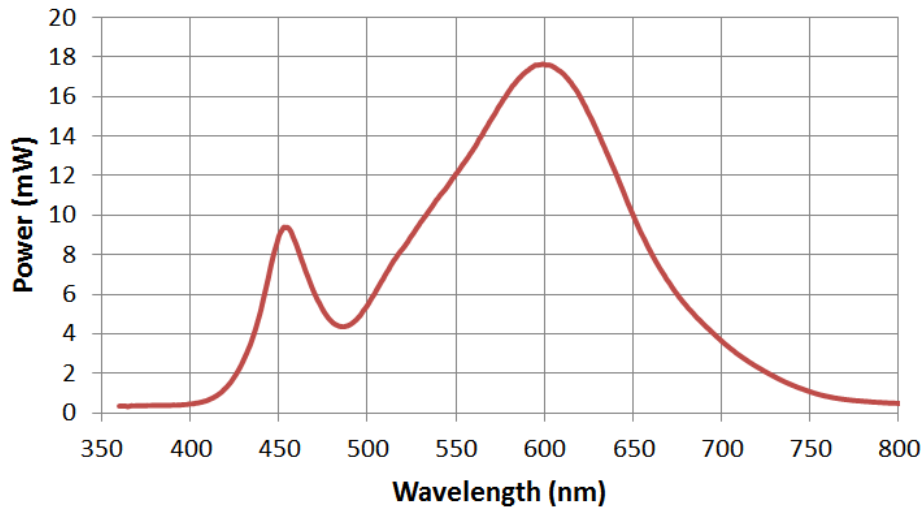
Figure 33: Typical relative spectral power vs wavelength (3000K model)

A20980/RGBW

Typical Output: 130 lumens (Red), 250 lumens (Green), 70 lumens (Blue), 400 lumens (White).

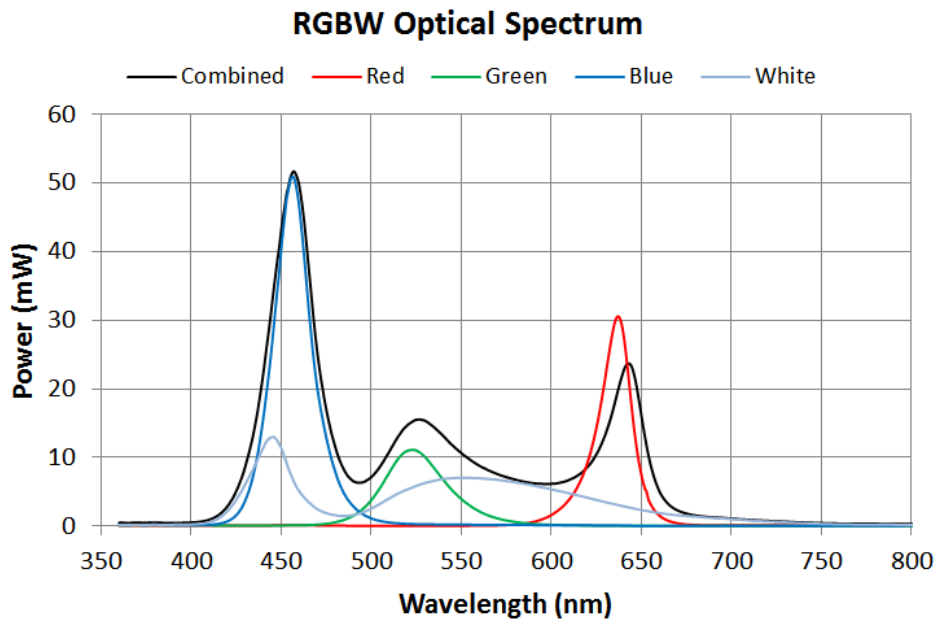
Figure 34: Typical relative spectral power vs wavelength (RGBW model)

Product Specification

SCHOTT
glass made of ideas

Replacement Parts

Multiport Terminal Blocks

Würth Elektronik (Part # **691 364 300 008**)

Power Supply

Mean Well USA (Part # **GS120A24-R7B**)

Fiber Optic Thumb Screw

McMaster Carr (Part # **94567A300**)

Rubber Feet

McMaster Carr (Part # **9541K8**)

Accessories

Locking USB Cable

Sentech USB (Part # **CBL-PD203MBS-2M**)

US Power Cable

TRC Electronics (Part # **3271V86**)

EU Power Cable

TRC Electronics (Part # **2171H576**)

UK Power Cable

TRC Electronics (Part # **210A71H576**)

Return Material Authorization (RMA)

Return Material Authorization (RMA)

Troubleshooting

If you are unsuccessful at resolving the following conditions, contact SCHOTT North America or your sales representative.

CAUTION

Do not attempt to repair the Light Source. Removing the cover will void the warranty.

Problem	Reason/Measure
Indicator LED on front panel does not light up.	Check for power connected to the Light Source.
Indicator LED on front panel is ON continuously, but LED output is dark.	Check polarity settings in software.
Indicator LED blinks	Fault exists. Use software interface to query reason.
Communications over Ethernet not possible even though network jack lights are flashing.	Ensure the Ethernet communication settings are correct.

Technical Support

If the Troubleshooting section does not resolve your issue, contact technical support at:
lightingimaging@us.schott.com

If technical support cannot help you resolve your issue remotely, you may need to return your unit under an RMA number. Technical support may suggest you elevate your RMA request to the engineering level. Make sure to mark this on your form to receive this service (additional fees may apply).

Return Material Authorization (RMA)

SCHOTT
glass made of ideas

Service and RMA Policy

Utilize the form on page 93 to gather information that will be needed by customer support. Contact customer support to receive an RMA number. Every field must be complete to receive more than a basic functionality repair and check. If any field is left blank, is not descriptive, or a problem is not reproducible, the repair technician will only repair the unit to pass our factory functional tester. If your problem is intermittent, you need to provide enough detail that the repair technician can replicate the issue to be able to repair it.

Any service required for any reason must be performed by SCHOTT or a SCHOTT authorized service representative. All services outside the warranty will be performed at the purchaser's request, according to normal service charges in effect at the time.

To return any item, whether for warranty repair or chargeable servicing, your RMA number must be clearly visible on the shipping label. All shipping charges must be prepaid.

SCHOTT guarantees all warranty repairs will be completed within two weeks of receipt. All units under warrantee will ship prepaid using our shipping method of choice. Alternate shipping methods will be shipped freight collect.

Warranty and Liabilities

SCHOTT warrants our light sources to be free from defective workmanship and materials. If, within two (2) years from shipment date, any product and/or part thereof are determined by SCHOTT to be defective, we will repair or replace it with new or reconditioned product and/or part.

Warranty is void if:

- We determine the product has been subjected to neglect or misuse, or has been installed following procedures not in accordance with our instruction manual.
- Unauthorized repairs or modifications have occurred.
- The warranty seal has been broken or the serial number label has been altered.

Our obligation is limited to repair or replacement. We will not be held responsible for consequential damages, transportation, installation, adjustment or other expenses arising in connection with our products or parts. This warranty is in lieu of all other statements or guarantees, written or implied, by SCHOTT or SCHOTT authorized representatives.

Return Material Authorization (RMA)

RMA Request Form

Note: All fields must be complete to receive an RMA number from customer service. When describing your problem, please include any information on which interface you are experiencing problems with, what you have tried to fix the issue, how to replicate the problem, etc.

Customer Information	
Name	
Company Name	
Return Shipping Address	
Phone	
Email	

CV-LS Information	
Serial Number	
Describe your complete setup and problem. Please include how the unit is powered, what controls you are using, ambient temperature, etc. Note: "Doesn't work" is not sufficient to diagnose your issue!	

Contact Information

SCHOTT
glass made of ideas

Customer Support

Lighting and Imaging
SCHOTT North America, Inc.
122 Charlton Street
Southbridge, MA 01550
Phone: (508) 765-9744
Fax: (508) 765-1299
E-Mail: lightingimaging@us.schott.com
Website: www.us.schott.com

Downloads

The data sheet, manuals, drivers, and the CV-LS Dashboard demo application can be downloaded from the SCHOTT website at the following link:

<https://www.us.schott.com/lightingimaging/english/sensors/products/coldvision/cvls.html>

Technical Support

You can reach us by e-mail at lightingimaging@us.schott.com

MicroscopeWorld

info@microscopeworld.com | 800-942-0528