# Instruction Manual ZEISS Primostar 1

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Manual de instrucciones ZEISS Primostar 1

## Instruction Manual ZEISS Primostar 1

Original Manual (en-US)

Translation of the original manual (zh-CHS, de, fr, it, pl, pt, ru, es)

Order Number: 415501-7021-111

Revision: 2

Language: en-US, zh-CHS, de, fr, it, pl, pt, ru, es

Effective from: 06/2021

Carl Zeiss Microscopy GmbH

Carl-Zeiss-Promenade 10

07745 Jena

Germany

[info.microscopy.de@zeiss.com](mailto:info.microscopy.de@zeiss.com)

[www.zeiss.com/microscopy](http://www.zeiss.com/microscopy)

Carl Zeiss Suzhou Co., Ltd.

Modern Industrial Square 3-B, No.333 XingPu Road SIP

215126 Suzhou

China

© 2021 Without the prior written consent of ZEISS, this document or any part of it must neither be translated nor reproduced or transmitted in any form or by any means - including electronic or mechanic methods, by photocopying, recording or by any information or filing system. The right to make backup-copies for archiving purposes shall remain unaffected thereby. Any violations may be prosecuted as copyright infringements.

The use of general descriptive names, registered names, trademarks, etc. in this document does not imply that such names are exempt from the relevant intellectual property laws and regulations and therefore free for general use. This shall also apply if this is not specifically referred to. Software programs shall entirely remain the property of ZEISS. No program or subsequent upgrade thereof may be disclosed to any third party, copied or reproduced in any other form without the prior written consent of ZEISS, even if these copies or reproductions are destined for internal use at the customer's only, the only exception being one single back-up copy for archiving purposes.

ZEISS reserves the right to make modifications to this document without notice.

# Table of Contents

|          |  |           |
|----------|--|-----------|
| <b>1</b> | <b>General Information.....</b>                                  | <b>5</b>  |
| 1.1      | Text Conventions and Link Types .....                            | 5         |
| 1.2      | Further Applicable Documents .....                               | 5         |
| 1.3      | Explanation of Warning messages and additional Information ..... | 6         |
| 1.4      | Contact.....   | 7         |
| <b>2</b> | <b>Safety .....</b>  | <b>8</b>  |
| 2.1      | Intended Use .....   | 8         |
| 2.2      | General Safety Information .....                                 | 8         |
| 2.2.1    | Requirements for Operators .....                                 | 8         |
| 2.2.2    | Safe Operating Condition.....                                    | 9         |
| 2.3      | Prevention of Hazards .....                                      | 9         |
| 2.3.1    | Mechanical Hazards .....   | 9         |
| 2.3.2    | Electrical Hazards .....   | 9         |
| 2.3.3    | Hazards Generated with the Operating Environment.....            | 9         |
| 2.3.4    | Ergonomic Hazards .....  | 10        |
| 2.3.5    | Hazards generated by Materials and Substances .....              | 10        |
| 2.4      | Warning Labels and Lights .....                                  | 10        |
| 2.4.1    | Meaning of Warning and Information Labels .....                  | 10        |
| 2.4.2    | Warning and Information Labels on the Stand .....                | 11        |
| <b>3</b> | <b>Product and Functional Description .....</b>                  | <b>12</b> |
| 3.1      | Main Components of Primostar 1.....                              | 12        |
| 3.2      | Control Elements and Indicators on the Stand.....                | 13        |
| 3.3      | Objectives .....   | 15        |
| 3.4      | Eyepiece.....  | 16        |
| <b>4</b> | <b>Installation.....</b>   | <b>17</b> |
| 4.1      | Connecting the Microscope to the Mains Power Supply.....         | 17        |
| <b>5</b> | <b>First Operating Steps .....</b>                               | <b>18</b> |
| 5.1      | Prerequisites for Commissioning and Operation .....              | 18        |
| 5.2      | Switching On the Microscope .....                                | 18        |
| 5.3      | Adjusting the Position of the Eyepieces .....                    | 18        |
| 5.4      | Installing the Blue Filter .....                                 | 19        |
| 5.5      | Acquiring an Image.....  | 19        |
| 5.6      | Changing Objectives .....  | 21        |
| 5.7      | Switching Off the Microscope .....                               | 21        |
| <b>6</b> | <b>Care and Cleaning Work .....</b>                              | <b>22</b> |
| 6.1      | Cleaning an Optical Surface .....                                | 23        |
| 6.2      | Removing Water-soluble Contamination .....                       | 23        |
| <b>7</b> | <b>Troubleshooting .....</b>                                     | <b>24</b> |

---

|           |  |           |
|-----------|--|-----------|
| <b>8</b>  | <b>Decommissioning and Disposal.....</b>                             | <b>25</b> |
| 8.1       | Decommissioning.....   | 25        |
| 8.2       | Decontamination .....  | 25        |
| 8.3       | Transport and Storage .....  | 25        |
| 8.4       | Disposal .....   | 26        |
| <b>9</b>  | <b>Technical Data and Conformity.....</b>                            | <b>27</b> |
| 9.1       | Applicable Standards and Regulations.....                            | 27        |
| 9.2       | Performance Data and Specifications .....                            | 27        |
| 9.3       | Performance Data and Specifications of the Optional Components ..... | 28        |
| <b>10</b> | <b>Accessories and Optional System Expansions .....</b>              | <b>30</b> |
| 10.1      | Installing the Mirror .....  | 30        |
|           | <b>Glossary .....</b>  | <b>31</b> |
|           | <b>Index.....</b>  | <b>32</b> |

# 1 General Information

This Instruction Manual is considered to be part of the Primostar 1. Herein after referred to as Microscope.

This Instruction Manual contains basic steps and safety information that must be observed during operation and maintenance. Therefore, the Instruction Manual must be read by the operator prior to commissioning and must always be available at the place of use of the Microscope. This Instruction Manual is an essential part of the Microscope and, if the Microscope is resold, the Instruction Manual must remain with the Microscope or be handed over to the new owner.

## 1.1 Text Conventions and Link Types

The following text conventions and link types are used:

| Text convention   | Meaning   |
|---|---|
| Click <b>Start</b> .<br>Press the <b>STANDBY</b> button.<br>Press <b>[Enter]</b> on the keyboard. | The names of controls and important information are shown in bold letters.  |
| Press <b>&lt;Ctrl+Alt+Del&gt;</b>   | Press several keys on the keyboard simultaneously.  |
| Select <b>Tools &gt; Goto Control Panel &gt; Airlock</b> .  | Follow a path in the software.  |
| <i>Text input</i>   | Text to be entered by the user  |
| Programming and Macros  | Anything typed in literally during programming, including, for example, macro codes, keywords, data types, method names, variables, class names, and interface names. |

Tab. 1: Text convention

| Link type   | Meaning                                     |
|---|---|
| See: <i>Text Conventions and Link Types</i> [▶ 5].  | Link to further information for this topic. |
| <a href="https://www.zeiss.com/corporate/int/home.html">https://www.zeiss.com/corporate/int/home.html</a> | Link to a website on the internet.          |

Tab. 2: Link types

## 1.2 Further Applicable Documents

Please take also note of the following documents:

|   |  |
|---|--|
| <b>Brochures and Certificates</b>                       | For Brochures, ISO certificates, CSA certificates and EU declarations of conformity ask your ZEISS Sales & Service Partner.  |
| <b>Local and National health and safety regulations</b> | Local and national health and safety regulations should be adhered to for the site and the use of the Microscope. Please consult with your ZEISS Sales & Service Partner in the event of these regulations conflicting with the installation requirements of the Microscope. |
| <b>Safety data sheets</b>                               | Observe the enclosed safety data sheets. The instructions and guidelines given in the respective safety data sheets must be complied with.   |
| <b>System and 3rd Party Components, Accessories</b>     | The Microscope can be configured in many ways. Information about the individual components, enhancements and accessories can be obtained from your ZEISS Sales & Service Partner. Also refer to the 3rd Party documentation of the manufacturer.                             |

## 1.3 Explanation of Warning messages and additional Information

DANGER, WARNING, CAUTION, and NOTICE are standard signal words used to determine the levels of hazards and risks of personal injury and property damage. Not only the safety and warning messages in the **Safety** chapter are to be considered also all safety and warning messages in other chapters. Failure to comply with these instructions and warnings can result in both personal injury and property damage and involve the loss of any claims for damages.

The following warning messages indicating dangerous situations and hazards are used in this document.

### **DANGER**

#### **Type and source of danger**

DANGER indicates an imminently hazardous situation which, if not avoided, will result in death or serious injury.

### **WARNING**

#### **Type and source of danger**

WARNING indicates a potentially hazardous situation which, if not avoided, may result in death or serious injury.

### **CAUTION**

#### **Type and source of danger**

CAUTION indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury.

### **NOTICE**

#### **Type and source of danger**

NOTICE indicates a potentially harmful situation which, if not avoided, may result in property damage.

### **Info**

Provides additional information or explanations to help users better understand the contents of this Instruction Manual.

## 1.4 Contact

If you have any questions or problems, please contact your local ZEISS Sales & Service Partner or one of the following addresses:

### Headquarters

|        |                              |
|--------|------------------------------|
| Phone: | +49 1803 33 63 34            |
| Fax:   | +49 3641 64 3439             |
| Email: | info.microscopy.de@zeiss.com |

### Courses and training

|        |                                 |
|--------|---------------------------------|
| Email: | courses.microscopy.de@zeiss.com |
|--------|---------------------------------|

### ZEISS Portal

The ZEISS Portal (<https://portal.zeiss.com/>) offers various services that simplify the daily work with your ZEISS systems (machines and software). It is constantly improved and extended to meet your needs and requirements better.

### ZEISS Sales & Service Partner

You can find a ZEISS Sales & Service Partner in your area under <https://www.zeiss.de/mikroskopie/website/forms/sales-and-service-contacts.html>.

### Service Germany

|        |                                 |
|--------|---------------------------------|
| Phone: | +49 7364 20 3800                |
| Fax:   | +49 7364 20 3226                |
| Email: | service.microscopy.de@zeiss.com |

## 2 Safety

This chapter contains general requirements for safe working practices. Any person using the Microscope or commissioned with installation or maintenance must read and observe these general safety instructions. Knowledge of basic safety instructions and requirements is a precondition for safe and fault-free operation. Operational safety of the supplied Microscope is only ensured if it is operated according to its intended use.

If any work is associated with residual risks, this is mentioned in the relevant parts of this document in a specific note. When components must be handled with special caution, they are marked with a warning label. These warnings must always be observed.

### 2.1 Intended Use

The Primostar 1 is a microscope that has been especially designed for education and routine.

Typical applications include:

- education (teaching, study)
- laboratories
- medical practices

Improper use can easily lead to impairment of its function or even damage to the Microscope. Damage caused by incorrect operation, negligence or unauthorized intervention, in particular by removing, modifying or replacing parts of the Microscope, can not be held liable by the manufacturer. Third-party devices or components that are not expressly approved by ZEISS may not be used.

### 2.2 General Safety Information

This Instruction Manual must be read before commissioning in order to ensure safe and uninterrupted operation. Pay particular attention to all listed safety notes. Make sure, that

- the operating personnel has read and understood this Instruction Manual, associated documents and particularly all safety regulations and instructions, and applies them.
- the local and national safety and accident prevention regulations must be observed, as well as the applicable laws and regulations in your country.
- this Instruction Manual is always available with/together with the Microscope.
- the Microscope is always in perfect condition.
- the Microscope is secured against access by unauthorized persons.
- maintenance and repair work, retrofitting, removal or replacement of components, as well as any other intervention in the Microscope not described in this Instruction Manual, may only be carried out by the manufacturer ZEISS or persons expressly authorized by ZEISS to do so.

#### 2.2.1 Requirements for Operators

The Microscope, components and accessories may only be operated and maintained by authorized and trained personnel. The Microscope may only be used in accordance this Instruction Manual. If the Microscope is not used as described, the safety of the user may be impaired and/or the Microscope may be damaged. Any unauthorized intervention or use other than within the scope of the intended use shall void all rights to warranty claims. The regional regulations on health protection and accident prevention must be observed at all times and during all work on and with the Microscope.

**Training** Authorized ZEISS personnel will provide basic training in operating the Microscope. As well as information on equipment safety and maintenance work that can be conducted by the operator. The training will be documented by ZEISS and its completion is to be confirmed by the operator.

Special application training is offered for a fee. Current training dates, additional information and the registration form can be found at <https://www.zeiss.com/microscopy/int/service-support/training-and-education.html>.

### 2.2.2 Safe Operating Condition

If circumstances occur which impair safety and cause changes in operating behavior, the Microscope must be shut down immediately and a ZEISS service representative should be informed. The Microscope can only be operated after proper installation by a ZEISS service representative and in compliance with the operating conditions. The Microscope may only be operated after correct installation by a ZEISS service representative and if the operating conditions are adhered to.

- Do not operate the Microscope until you have completely read and understood the entire documentation delivered with the Microscope.
- Make sure that all protective cover panels are installed and all warning labels are available and legible.
- Ensure conditions and take measures to prevent the build up of electrostatic charge on the workplace.

## 2.3 Prevention of Hazards

This section summarizes potential hazards and recommended safety precautions. Failure to follow the safety instructions and instructions may result in personal injury and property damage.

### 2.3.1 Mechanical Hazards

**Property Damage due to Transport** There is a risk of injury and property damage if the Microscope is improperly handled and transported.

- Only use the handle, if applicable, for transport of the Microscope. Otherwise hold the Microscope with one hand and the base plate with the other hand.

### 2.3.2 Electrical Hazards

**Voltage Hazards** Risk of electric shock in case of contact with live parts.

Always use the power adapter supplied by ZEISS. When an unsuitable power adapter is used, ZEISS can no longer guarantee the electrical safety and functionality of the Microscope.

- Shut down the Microscope.
- Disconnect the power supply before cleaning.
- Set up and operate the Microscope so that the connectors are easily accessible.

### 2.3.3 Hazards Generated with the Operating Environment

**Dirt, Dust and Moisture** Dirt, dust and moisture can impair the Microscope's functionality.

- Shut down the Microscope whenever it is not used and cover it with a dust protection cover.
- Always cover unused openings/ports.
- Regular maintenance and cleaning according to the instructions in this document.
- Make sure that no cleaning liquid or moisture gets inside the Microscope.
- Make sure that the electrical parts never come into contact with moisture.
- Never expose the Microscope to inadmissible climate conditions (high humidity and temperature).

### 2.3.4 Ergonomic Hazards

**Prevention of Musculoskeletal Disorders** Musculoskeletal disorders (MSDs) affect the muscles, nerves, blood vessels, ligaments and tendons. Workers in many different industries and occupations can be exposed to risk factors at work, such as lifting heavy items, bending, reaching overhead, pushing and pulling heavy loads, working in awkward body postures and performing the same or similar tasks repetitively. Employers are responsible for providing a safe and healthful workplace for their workers.

### 2.3.5 Hazards generated by Materials and Substances

- Infection Hazards** Direct contact with the eyepieces can be a potential way of passing on bacterial and viral infections.
- The risk can be lowered by using personal eyepieces or eyecups. If eyepieces need to be disinfected frequently, ZEISS recommends to use the eyepieces without eyecups.
  - To avoid infections, the use of personal protective equipment (PPE) e.g. gloves for operation, cleaning, and decontamination is highly recommended. Disposable gloves can be decontaminated with, e.g., alcohol, if necessary, or should be changed frequently to minimize the risk of contamination.
- Consumable Hazards** Incorrect handling of consumables and cleaning agents can lead to property damage or skin and eye injuries. Consumables that are not approved by ZEISS can lead to property damage. Consult your ZEISS Sales & Service Partner to learn what consumables you can order and how to handle them.
- Disinfectant Hazards** Ensure adequate ventilation in closed rooms. In case of insufficient ventilation, wear respiratory protective equipment. Remove any harmful residue. Allow the device to dry off after disinfection, particularly after disinfection of eyepieces. Do not inhale vapors. When using disinfectants, do not eat, drink or smoke. Avoid contact with eyes and skin. Remove contaminated clothing and wash before reuse.
- Eye, Skin, Respiratory Tract Irritation** Exposure to chemicals and their aerosols can cause eye, skin and respiratory tract irritation. Use appropriate personal protective equipment (PPE).

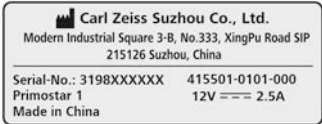
## 2.4 Warning Labels and Lights

All parts of the Microscope that may pose specific hazards are marked with additional warning labels (pictograms) on the Microscope. These warning labels indicate potential hazards and form part of this Instruction Manual. They are to be kept in clean and easily legible condition.

Check all the mandatory warning labels for: Availability, Legibility, Correctness. Damaged or illegible warning labels must be replaced immediately. Always observe all warning labels on the complete Microscope.

### 2.4.1 Meaning of Warning and Information Labels

The meaning of the individual warning labels is explained in the following:

| Pos. | Symbol  | Description           |
|------|---|-----------------------|
| 1    |  <p>Carl Zeiss Suzhou Co., Ltd.<br/>Modern Industrial Square 3-B, No.333, XingPu Road SIP<br/>215126 Suzhou, China<br/>Serial-No.: 3198XXXXXX 415501-0101-000<br/>Primostar 1 12V --- 2.5A<br/>Made in China</p> | Microscope type label |

| Pos. | Symbol  | Description   |
|------|---|---|
| 2    |  | Hot surface below<br>Operate microscope only on a stable, solid, smooth and not tinderlike (non-flammable) surface. |
| 3    |  | Microscope type label   |
| 4    |  | Primostar 1<br>EU Importer<br>Carl Zeiss Microscopy GmbH<br>Carl-Zeiss-Promenade 10<br>07745 Jena, Germany          |

2.4.2 Warning and Information Labels on the Stand

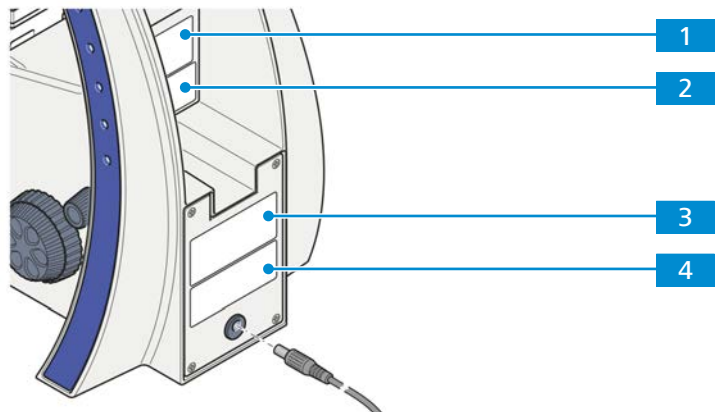
Fig. 1: Position of warning and information labels

## 3 Product and Functional Description

The Primostar 1 is a transmitted-light microscope of compact design with a small footprint.

The Primostar 1 is a microscope that has been specially developed for education and routine. It features great durability in permanent use.

### 3.1 Main Components of Primostar 1

Primostar 1 comprises the following main components

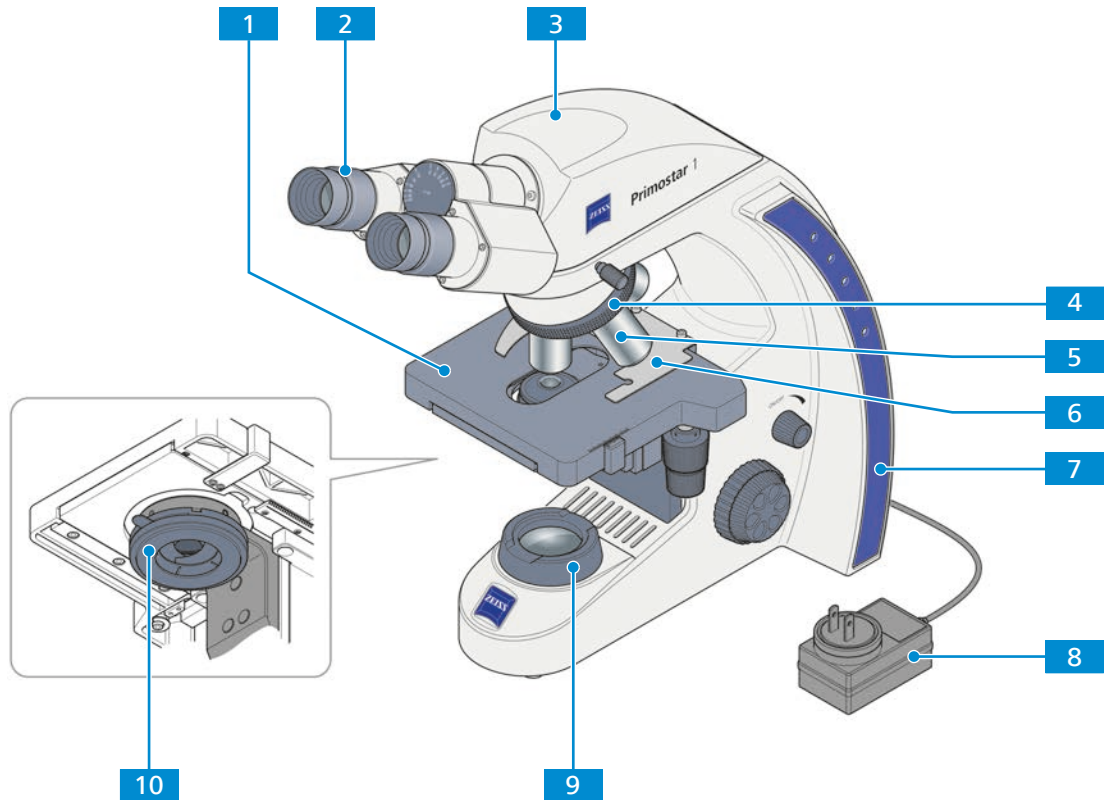
Fig. 2: Main Components of Primostar 1

- |  |  |
|--|--|
| <b>1</b> Microscope stage                    | <b>2</b> Eyepiece                      |
| <b>3</b> Tube                                | <b>4</b> Nosepiece                     |
| <b>5</b> Objective                           | <b>6</b> Sample holder                 |
| <b>7</b> Stand                               | <b>8</b> Power supply unit             |
| <b>9</b> Transmitted-light illumination unit | <b>10</b> Abbe condenser, Fixed Köhler |

### 3.2 Control Elements and Indicators on the Stand

**Purpose** The controls on the stand control the main functions of the microscope.

**Position** The following controls are located on the right side of the stand.

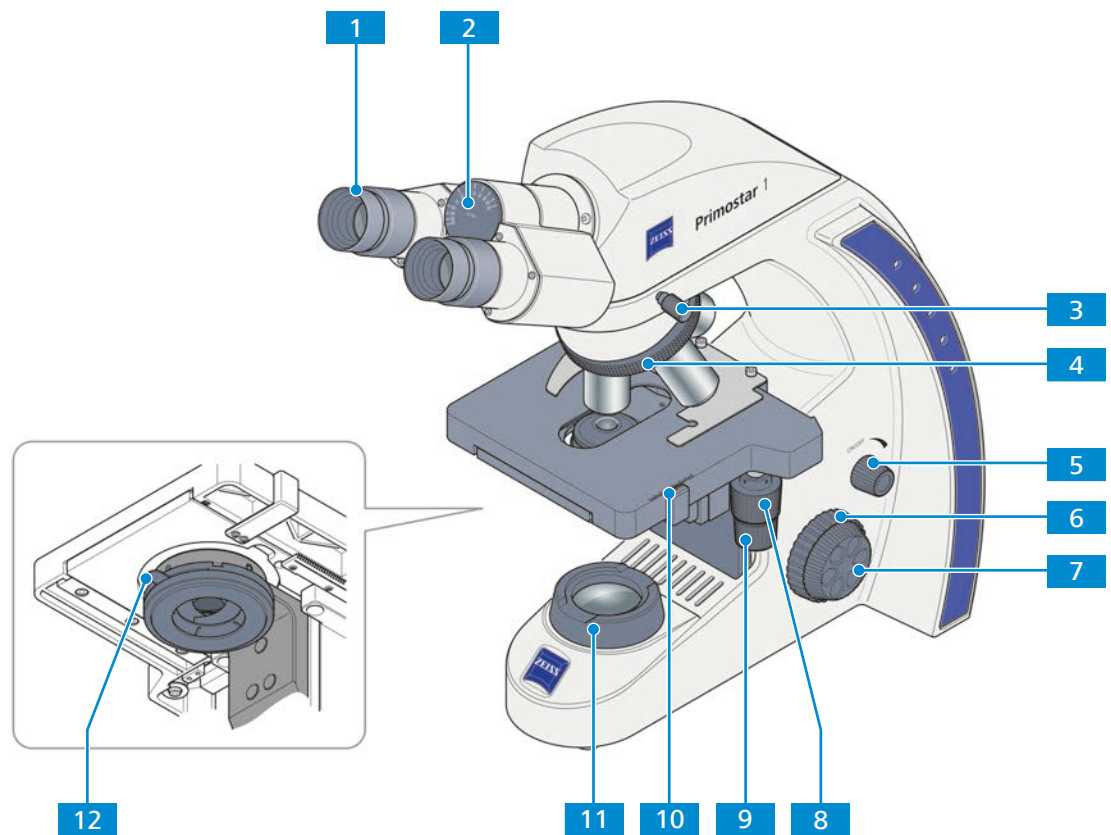
Fig. 3: Stand controls on the right side

- |  |   |
|--|---|
| <b>1</b> Eyepiece, adjustable                        | <b>2</b> Interpupillary distance indicator                            |
| <b>3</b> Fixing screw observation tube               | <b>4</b> Knurled ring for turning the nosepiece                       |
| <b>5</b> Light intensity knob                        | <b>6</b> Coarse focusing drive (right side)                           |
| <b>7</b> Fine focusing drive (right side)            | <b>8</b> Control knob for Y travel of rackless stage                  |
| <b>9</b> Control knob for X travel of rackless stage | <b>10</b> Vernier and scale, displaying the Y position of the stage   |
| <b>11</b> Transmitted-light illumination unit        | <b>12</b> Lever for adjusting the aperture diaphragm of the condenser |

**Position** The following controls are located on the left side of the stand.


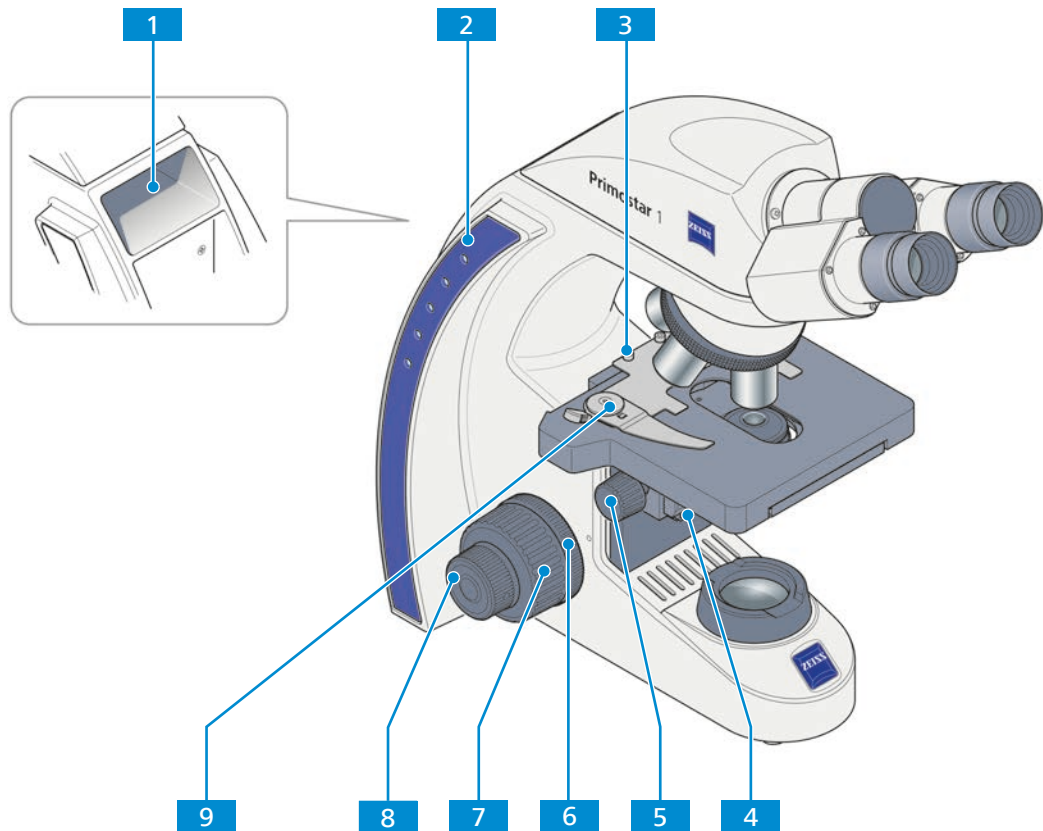
Fig. 4: Stand controls on the left side

- |   |   |
|---|---|
| <b>1</b> Carrying handle                                  | <b>2</b> Illumination-intensity indicators for transmitted light                |
| <b>3</b> Clamping screw for the sample holder             | <b>4</b> Fixing screw for lowering the condenser                                |
| <b>5</b> Knurled ring for condenser height adjustment     | <b>6</b> Knurled ring for adjusting the smoothness of the coarse focusing drive |
| <b>7</b> Coarse focusing drive (left side)                | <b>8</b> Fine focusing drive (left side)  |
| <b>9</b> Lever of the sample holder for fixing the sample |   |

### 3.3 Objectives

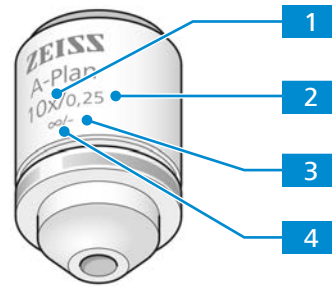The functional characteristics of an objective are indicated by the respective labeling, e. g. **A-Plan 10x/0,25 ∞/-**.

#### 1 Magnification of the objective

The magnification factor specifies the magnification of the objective. Larger magnifications allow to resolve smaller features on the sample.

The objective's magnification factor is also expressed by the objective's color ring, see ZEISS color code.

The objective magnification multiplied by the eyepiece magnification gives the overall magnification.

#### 2 Numerical aperture (NA)

The numerical aperture multiplied by 1000 represents the maximum useful magnification - above that limit no additional resolution is provided.

#### 3 Applicable cover glass thickness

- : The objective can be used without cover glass or with a cover glass with a thickness of 0.17 mm.

0: The objective can be used without cover glass.

0.17: The objective can be used with a cover glass with a thickness 0.17 mm.

#### 4 Mechanical tube length

The objective can be used with tubes of any length.

#### Info

Due to their short working distance, the following objectives have a resilient mount (sample protection):

- ▶ 40x
- ▶ 100x oil

### 3.4 Eyepiece

**Purpose** The eyepieces serve to observe the microscopic image.

**Position** The eyepieces are inserted into the tube.

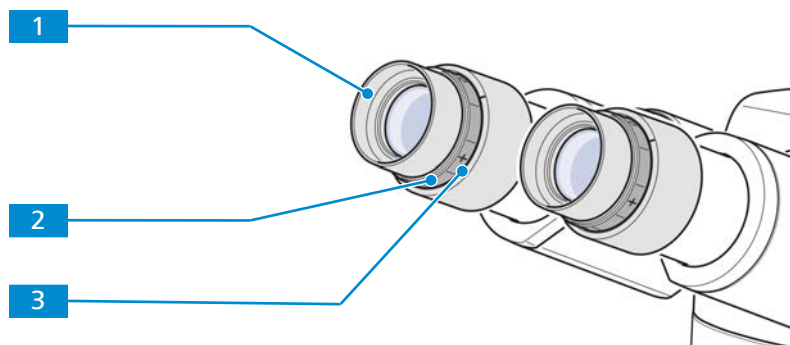
Fig. 5: Eyepiece

**1** Foldover rubber eyecup

**2** Focusing ring  
for compensation of defective vision

**3** Diopter scale  
to facilitate finding the correct setting

**Function** Both eyepieces are suitable for spectacle wearers. Additionally, they contain a focusing ring for compensation of defective vision. The provided diopter scale helps to find the correct setting.


## 4 Installation

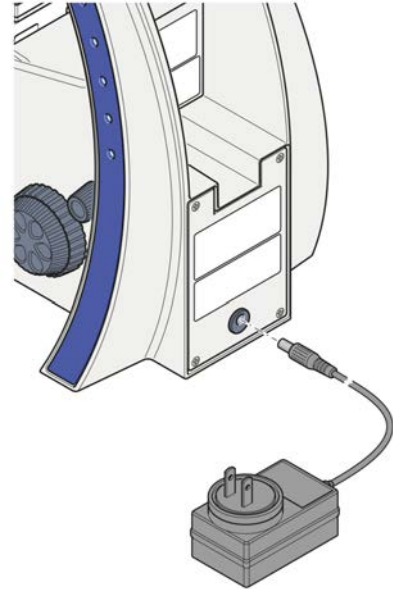Perform only the installation work described in this document. All other installation work not described below may only be carried out by an authorized ZEISS service representative.

### 4.1 Connecting the Microscope to the Mains Power Supply

- Procedure**
1. Remove the plug-in power unit from its storage fixture.
  2. Connect the plug-in power unit to the mains socket of the microscope.

3. If necessary, replace the installed power outlet adapter with one of the supplied country-specific adapters.
4. Connect the plug-in power unit to the mains power supply.

## 5 First Operating Steps

This chapter describes the switching on/off as well as the first operating steps with the Microscope.

### Info

Additional information and detailed descriptions are available in the further applicable documents, or ask your ZEISS Sales & Service Partner.

### Info

Further information on the software and its operation is available in the software's online help.

### 5.1 Prerequisites for Commissioning and Operation

The following basic prerequisites are necessary for commissioning and operation:

- The Instruction Manual was read prior to commissioning or operation and kept for further use.
- Chapter **Safety** is read and understood.
- Operator is acquainted with the general Windows®-based programs.
- If required: Basic training and safety briefing successfully completed.

### 5.2 Switching On the Microscope

**Prerequisite** ✓ *The microscope is connected to the mains [▶ 17].*

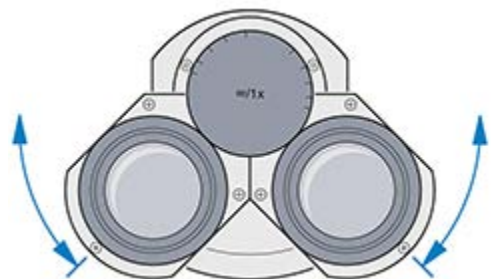
- Procedure**
1. Turn the light intensity knob.
  2. Adjust the illumination to the desired intensity.
    - The selected intensity is indicated by the light-emitting diodes on the stand.

### 5.3 Adjusting the Position of the Eyepieces


### Info

The adjustment of the interpupillary distance is correct when you see only one round image while looking through the two eyepieces.

- Procedure**
1. Set the interpupillary distance by rotating the eyepiece tubes symmetrically toward or away from one another.

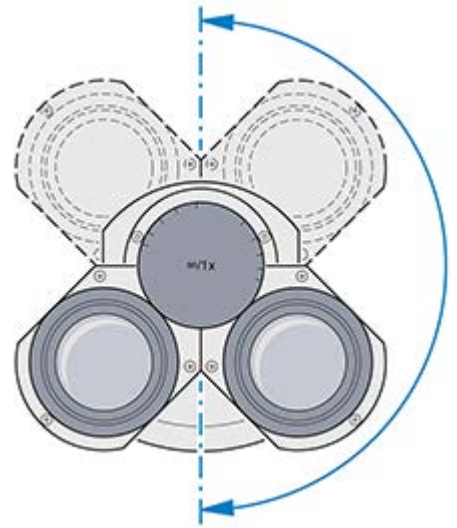
2. Set the viewing height by swivelling the eyepieces a full 180° upwards or downwards.

## 5.4 Installing the Blue Filter

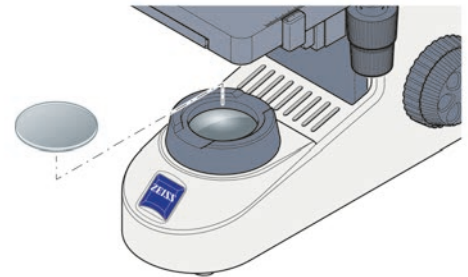- Procedure**
1. Put the blue filter onto the luminous-field diaphragm.

Proceed in the reverse order for removal.


## 5.5 Acquiring an Image
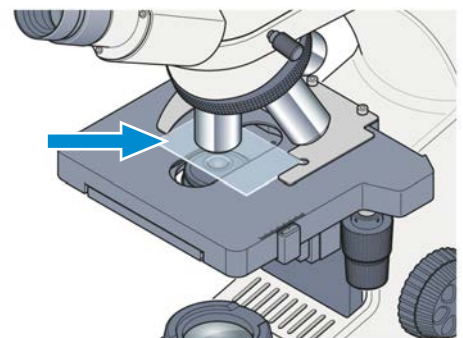
### Info


The microscope is supplied factory-adjusted

**Prerequisite** ✓ *The position of the eyepieces is set [▶ 18].*


- Procedure**
1. Place the sample in the sample holder of the mechanical stage.

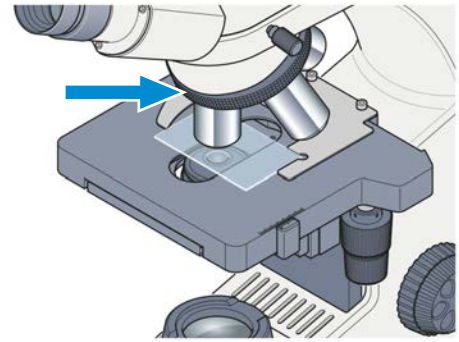
2. Select the desired magnification by placing the corresponding objective in the light path.

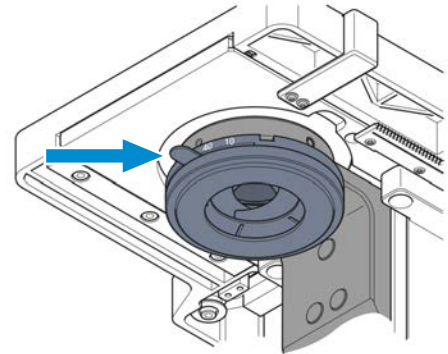
3. Set the control lever of the condenser aperture diaphragm to the value of the selected magnification (**10x**, **40x** or **100x**).

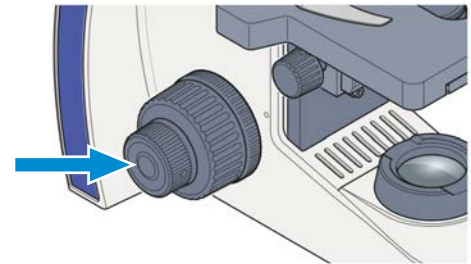4. Focus on the sample using the focusing drive.

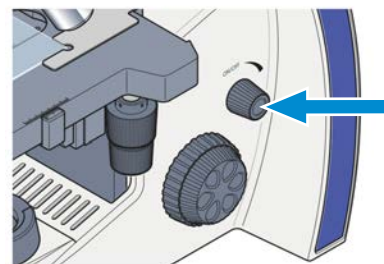5. Use the rotary knob for illumination intensity to adjust the illumination to a comfortable setting.

## 5.6 Changing Objectives

### NOTICE

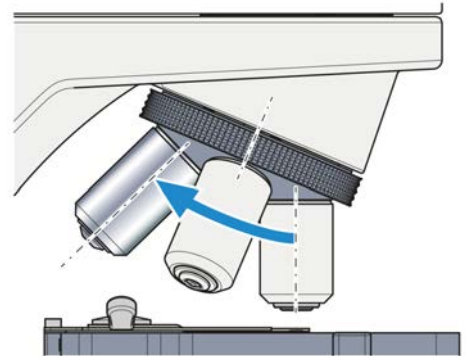
#### Dust-sensitive components

If unused nosepiece openings remain uncovered, particles may enter the Microscope and may damage its optics and mechanics permanently.


- ▶ Always close unused nosepiece openings with cover caps!

The objectives should be installed according to the direction shown in the figure in order of increasing magnification.

- Procedure**
1. Use the focus drive to fully lower the mechanical stage or the stage carrier.
  2. Remove the cover cap or screw out the present objective from the nosepiece's relevant opening.

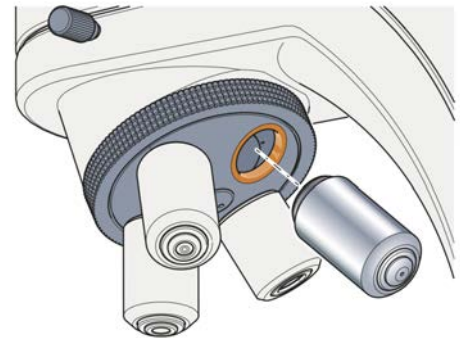3. Take the objective out of its case.
4. Carefully screw the objective into the opening. Make sure it engages properly in the nosepiece's thread.

## 5.7 Switching Off the Microscope

- Procedure**
1. Turn the light intensity knob.
    - The light-emitting diodes turn off.
  2. Cover the microscope with the dust cover.

## 6 Care and Cleaning Work

To ensure the best possible performance of the Microscope, maintenance must be performed on a regular basis. To maintain operational safety and reliability of the Microscope, we recommend entering into a **ZEISS Protect service agreement**. Please keep the service logs for your Microscope.

### **DANGER**

#### **Electric shock due to live parts**

When the Microscope is still switched on, coming in contact with live parts can lead to electric shock or burn.

- ▶ Switch off Microscope prior to opening or cleaning.
- ▶ Disconnect live parts from the power supply.

### **NOTICE**

#### **Functional impairment due to dirt and moisture**

Dirt, dust and moisture can impair the Microscope's functionality and can cause short-circuits.

- ▶ Use the dust protection cover if the Microscope is not used.
- ▶ The ventilation slots must be unobstructed at all times.
- ▶ Perform regular maintenance and cleaning according to the instructions in this document and according to the instructions in the applicable documents.
- ▶ Make sure that no cleaning liquid or moisture gets inside the Microscope.
- ▶ In case of damage, the affected parts of the Microscope must be taken out of operation.

## 6.1 Cleaning an Optical Surface

### NOTICE

#### Damage of the lens due to improper cleaning

- ▶ Remove dust from the lens slowly and carefully.
- ▶ Remove dust on optical surfaces with a natural-hair brush or blow it off with a rubber bellows.

- Parts and Tools**
- 🔧 Clean cloth
  - 🔧 Cotton swab
  - 🔧 Optical cleaning solution (85% n-hexan and 15 vol% isopropyl alcohol (IPA))
  - 🔧 Lint-free cloth

- Procedure**
1. Moisten a cotton swab or a clean cloth with an optical cleaning solution, if necessary.
  2. Wipe optical surfaces in a circular motion towards the edge of the optics with slight pressure.

WRONG

CORRECT

3. Dry with a lint-free cloth.

## 6.2 Removing Water-soluble Contamination

- Parts and Tools**
- 🔧 Clean cloth
  - 🔧 Lint-free cloth

- Procedure**
1. Moisten a clean cloth with water.  
→ A mild detergent may be added to the water (no solvent!).
  2. Wipe off the area with the cloth.
  3. Dry with a lint-free cloth.

## 7 Troubleshooting

The following table provides information about solving common problems.

### Info

If you cannot solve the problem or if you are unsure about a certain technical difficulty, contact your local ZEISS service representative.

| Symptom  | Cause   | Measure   |
|--|---|---|
| The field of view is not completely visible                                | Nosepiece with objective is not completely switched into the click-stop position.     | Switch the nosepiece with objective into the click-stop position.   |
|  | The filter is not placed correctly on the luminous-field diaphragm.                   | <i>Place the filter correctly</i> [▶ 19].   |
| Low resolving power and poor image contrast                                | The aperture diaphragm is not correctly adjusted.                                     | Set the aperture diaphragm according to the 2/3 rule or the sample features.  |
|  | The condenser is not correctly focused.   | Focus the condenser.  |
|  | A wrong cover glass thickness for transmitted light objectives is used.               | Use standard cover glass with a thickness of 0.17 mm if objectives are corrected for 0.17 mm cover glass thickness. |
|  | Immersion objectives are used with no or inappropriate immersion oil.                 | Use immersion oil 518 N or 518 F from ZEISS.  |
|  | The immersion oil contains air bubbles.   | Repeat the oiling procedure with fresh oil.   |
|  | The front lens of a dry objective is soiled with immersion oil.                       | Clean the front lens of the dry objective.  |
|  | Dirt or dust on the optical surfaces of objectives, eyepieces, condensers or filters. | Clean the respective optical components.  |
| Significant difference in the focus position after changing the objective. | The focusing eyepieces are not set correctly.   | Set the focusing eyepieces to the corresponding defective vision.   |
| The LED source does not light when turning the light intensity knob.       | The power plug is not plugged into the power outlet.                                  | <i>Connect the power plug to the power outlet</i> [▶ 17].   |
| The stage comes down by itself, the focus is unstable.                     | The adjusted torque of the coarse focusing drive is too low.                          | Contact the ZEISS service representative.   |

## 8 Decommissioning and Disposal

This chapter contains information on the decommissioning and disposal of the Microscope and its expansions/components or accessories.

### 8.1 Decommissioning

If the Microscope is not used for an extended period such as several months, it should be shut down completely and secured against unauthorized access. Complete decommissioning of the Microscope should be executed by your ZEISS service representative.

#### **DANGER**

##### **Electric shock due to live parts**

When the Microscope is still switched on, coming in contact with live parts can lead to electric shock or burn.

- ▶ Switch off Microscope prior to opening or cleaning.
- ▶ Disconnect live parts from the power supply.

- Procedure**
1. Switch off the Microscope.
  2. Pull the mains plug.

### 8.2 Decontamination

A decontamination statement must be submitted before returning any used objects to the ZEISS location.

If reliable decontamination cannot be guaranteed, the hazard must be marked according to applicable regulations. In general, a well-visible warning sign must be affixed to the article itself and to the outside of the packaging, together with detailed information on the type of contamination.

### 8.3 Transport and Storage

The following regulations must be observed before and during transport:

- The boxes must be secured during transport.
  - Avoid rocking the boxes back and forth.
  - Note the weight information on the package and on the shipping document.
  - Where possible, the original packaging must be used for shipping or transport.
- Maximum shock resistance**
- Do not drop or bump the boxes during movement or storage. Any acceleration shall be < 10 g.
  - Evaluate packaging shock and tilting sensors on delivery and after internal transport.
- Allowable temperature**
- Allowable temperature during on-site storage:
- Between -10 °C and +40 °C
  - Relative humidity less than 80 % at +40 °C
- Allowable temperature during transport in packaging:
- Between -40 °C and +70 °C

**Info**

**24 hours before installation** of the Microscope it is required that the boxes be at recommended room temperature to avoid ingress of humidity, which is very harmful to optical paths, and to ensure effective stability of the Microscope during installation and testing.

## 8.4 Disposal

The Microscope and its components must not be disposed of as domestic waste or through municipal disposal companies. They must be disposed of in accordance with applicable regulations (WEEE Directive 2012/19/EU). ZEISS has implemented a system for the return and recycling of devices in member states of the European Union that ensures suitable reuse according to the EU Directives mentioned. The customer is responsible for decontamination.

**Info**

Detailed information on disposal and recycling is available from your ZEISS Sales & Service Partner.

## 9 Technical Data and Conformity

This chapter contains important technical data as well as information on the conformity.

### 9.1 Applicable Standards and Regulations

Observe all general and country-specific safety regulations as well as applicable environmental protection laws and regulations.

The Microscope is in compliance with the requirements of the following regulations and directives:

|   |   |
|---|---|
| 2011/65/EU<br>2015/836/EU               | RoHS Directive  |
| 98/79/EC                                | IvD Directive   |
| EN 55011                                | Noise emission according to CISPR 11 Group 1 Class B intended use in industrial environment   |
| EN 61010-1 and IEC 61010-2-101          | Safety requirements for electrical equipment for measurement, control, and laboratory use   |
| DIN EN 61326-1<br>DIN EN 61326-2-6:2013 | Electrical equipment for measurement, control and laboratory use – EMC requirements / IVD requirements  |
| EN 62471                                | Photobiological safety of lamps and lamp systems<br>Compliance with the limit values in accordance with EN 62471 Risk group 1 (Exempt risk group viewing from ocular) |

The Microscope and its accessories have been classified as instrument category 9 (laboratory equipment or comparable standard). It also complies, as applicable, with the EU-regulations 2011/65/EU (RoHS) and 2012/19/EU (WEEE).

European and International Directives / Standards: For more information on ISO and CSA certificates or CE Declarations of Conformity, contact your ZEISS Sales & Service Partner.

ZEISS works according to a certified Environment Management System according to ISO 14001. The Microscope was developed, tested and produced in accordance with the valid regulations and guidelines for environmental law of the European Union.

### 9.2 Performance Data and Specifications

The Microscope must only be operated in closed rooms. It is recommended to install the Microscope in a dark room where artificial illumination, sunlight or other light sources cannot interfere with image acquisition. The Microscope should not be installed near windows with direct sunlight or radiators. Compliance with the installation requirements of the Microscope and the availability of the requested supplies is the responsibility of the customer and has to be provided at the time of installation. Due to continuous development, we reserve the right to change specifications without notice.

The Microscope must be plugged into a properly installed power socket with protective earth contact using the supplied mains cable. The protective earth connection must not be impaired by the use of extension cables.

#### Info

Your ZEISS Sales & Service Partner will provide you with the detailed installation requirements.

| <b>Weight and Sizes</b>             | <b>Main Components</b>          | <b>Length (mm)</b>                         | <b>Width (mm)</b> | <b>Height (mm)</b> | <b>Weight (kg)</b> |
|-------------------------------------|---------------------------------|--|-------------------|--------------------|--------------------|
|                                     | Stand with binocular tube       | 410  | 190               | 395                | 7.7                |
| <b>Air Conditioning and Quality</b> | Temperature range for operation | 5 to 40 °C                                 |                   |                    |                    |
|                                     | Relative humidity               | < 80 % at 35 °C                            |                   |                    |                    |
|                                     | Atmospheric pressure / altitude | 800 to 1060 hPa / ≤ 2000 m above sea level |                   |                    |                    |
|                                     | Pollution degree                | 2  |                   |                    |                    |
| <b>Plug-in power unit</b>           | Nominal AC voltage              | L/N 100 to 240 V ± 10 %                    |                   |                    |                    |
|                                     | Nominal frequency               | 50 / 60 Hz                                 |                   |                    |                    |
|                                     | Main Power Plug                 | Local mains plug will be supplied.         |                   |                    |                    |
|                                     | Power consumption               | max. 0.8 A                                 |                   |                    |                    |
|                                     | Output                          | 12 V DC, max. 2.5 A                        |                   |                    |                    |
| <b>Mains connection</b>             | Stand rated input               | 12 V DC, 2.5 A                             |                   |                    |                    |
|                                     | Protection class                | IP20 (IEC 60529)                           |                   |                    |                    |
|                                     | IEC earth class                 | Class III of IEC 61140                     |                   |                    |                    |
|                                     | Overvoltage Category            | II   |                   |                    |                    |

### 9.3 Performance Data and Specifications of the Optional Components

The customer is responsible for ensuring that the installation conditions for the microscope are met and that the required equipment is already available at the time of installation. Changes are reserved due to continuous technical developments.

| <b>LED illumination</b>                            |  |                     |
|--|--|---------------------|
| LED  |  | white light         |
| Peak wavelength                                    |  | 440 nm              |
| LED class  |  | 1                   |
| Constant, brightness-independent color temperature |  | 3200 K              |
| Homogeneous field illumination                     |  | 20 mm (diameter)    |
| Suitable for objectives with magnifications of     |  | 4x to 100x          |
| Analogous brightness adjustment from               |  | approx. 15 to 100 % |

|                                  |   |  |
|----------------------------------|---|--|
| <b>Stand with stage focusing</b> | Coarse focusing drive                       | 42 mm / rev.                             |
|                                  | Fine focusing drive                         | 0.2 mm / rev.                            |
|                                  | Total stage lift                            | 15 mm                                    |
| <b>Nosepiece</b>                 | Objective change                            | manual via quadruple objective nosepiece |
|                                  | Objectives                                  | infinity-corrected objective range       |
|                                  | Mounting thread                             | W 0.8                                    |
| <b>Eyepieces</b>                 | Tube size                                   | 30 mm                                    |
|                                  | Field-of-view number                        | 20                                       |
|                                  | Magnification                               | 10x                                      |
|                                  | Suitable for spectacle wearers              | Br.                                      |
|                                  | Focusable                                   | Foc.                                     |
| <b>Sample stage</b>              | Type  | Rackless                                 |
|                                  | Dimensions (width x depth)                  | 140 x 140 mm                             |
|                                  | Stage travel (X x Y)                        | 75 x 40 mm                               |
|                                  | Coaxial drive                               | optional right or left                   |
|                                  | Vernier scales                              | readable from the right                  |
|                                  | Sample holder                               | with spring lever, left                  |
| <b>Condenser</b>                 | Abbe condenser 0.9; Fixed-Köhler            | for $V_{obj}$ 4x to 100x                 |
| <b>Illuminating mirror</b>       | Plane surface and spherical surface with f' | 75 mm                                    |
| <b>Binocular tube 30°/20</b>     | Maximum field-of-view number                | 20                                       |
|                                  | Interpupillary distance                     | adjustable from 50 to 75 mm              |
|                                  | Tube angle                                  | 30°                                      |
|                                  | Viewing height                              | 380 to 415 mm                            |
|                                  | Viewing port, tube factor                   | 1x                                       |

## 10 Accessories and Optional System Expansions

Only the following accessories may be used with the Microscope as their safe use has been confirmed by ZEISS. Only original parts from ZEISS may be used. Check in advance whether your Microscope can be retrofitted with a system expansion.

After installation or conversion it must be carefully checked whether the Microscope and its system expansions/components are in a safe operational state and whether unused ports on the stand are closed. For details and safety measures please refer to the associated documents.

### Info


Additional information and detailed descriptions are available in the further applicable documents, or ask your ZEISS Sales & Service Partner.

| Name       | Description/Info  |
|------------|---|
| Objectives | <p>The performance of the microscope objectives affects the image quality of your microscope like no other system component. Whether you work with histological samples, cell samples or entire organisms – the suitability of microscope objectives for your application depends on various factors.</p> <p>More detailed information on available and recommended objectives can be found at <a href="https://www.micro-shop.zeiss.com/de/de/shop/objectives">https://www.micro-shop.zeiss.com/de/de/shop/objectives</a> or ask your ZEISS Sales &amp; Service Partner.</p> |
| Mirror     | When no power outlet is available, the mirror serves to illuminate the sample.  |

### 10.1 Installing the Mirror

**Function** The mirror serves to illuminate the sample if no power outlet is available.

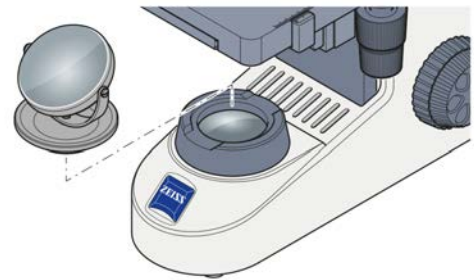**Procedure** 1. Place the mirror onto the mount of the luminous-field diaphragm.

2. Rotate and incline the mirror until the daylight is reflected homogeneously into the light path.

Proceed in the reverse order for removal.

# Glossary

**NA**

---

Numerical Aperture

**User**

---

Person examining a sample under the microscope.

**ZEISS**

---

ZEISS is an internationally leading technology enterprise operating in the fields of optics and optoelectronics. Further information about ZEISS can be found at [www.zeiss.com](http://www.zeiss.com).

**ZEISS Sales & Service Partner**

---

The Sales & Service Partner is generally in the field for customer support in a regional area and / or a clearly defined customer group.

**ZEISS service representative**

---

Specially trained service expert, either ZEISS staff or authorized service partner of ZEISS.

# Index

## A

|                              |    |
|------------------------------|----|
| Accessories                  | 30 |
| Air Conditioning and Quality | 28 |

## B

|             |    |
|-------------|----|
| Brightfield | 19 |
|-------------|----|

## C

|                              |    |
|------------------------------|----|
| Cleaning                     |    |
| Water-soluble contaminations | 23 |
| Connecting the microscope    | 17 |
| Contamination                | 25 |
| Controls                     | 13 |

## D

|                 |    |
|-----------------|----|
| Decontamination | 25 |
| Disposal        | 26 |

## E

|          |    |
|----------|----|
| Eyepiece | 16 |
|----------|----|

## F

|        |    |
|--------|----|
| Filter | 19 |
|--------|----|

## G

|                            |   |
|----------------------------|---|
| General Safety Information | 8 |
|----------------------------|---|

## H

|            |   |
|------------|---|
| Hazards    | 9 |
| Prevention | 9 |

## I

|                         |    |
|-------------------------|----|
| Image                   | 19 |
| Information labels      | 11 |
| Installing              |    |
| Filter                  | 19 |
| Mirror                  | 30 |
| Objective               | 21 |
| Intended Use            | 8  |
| Interpupillary distance | 18 |

## M

|                  |    |
|------------------|----|
| Mains connection | 28 |
| Maintenance      | 22 |
| Mirror           | 30 |

## N

|                    |    |
|--------------------|----|
| Numerical aperture | 15 |
|--------------------|----|

## O

|                            |        |
|----------------------------|--------|
| Objective                  | 15, 21 |
| Operation                  |        |
| Prerequisites              | 18     |
| Optional system expansions | 30     |
| Installation               | 30     |

## P

|                  |    |
|------------------|----|
| Performance data | 27 |
| Prerequisites    |    |
| Operation        | 18 |

## R

|               |   |
|---------------|---|
| Requirements  |   |
| for Operators | 8 |

## S

|                          |    |
|--------------------------|----|
| Safe Operating Condition | 9  |
| Safety                   | 8  |
| Shutdown                 | 25 |
| Switching off            | 21 |
| Switching on             | 18 |

## T

|                               |    |
|-------------------------------|----|
| Training                      | 8  |
| Transmitted-light brightfield | 19 |
| Troubleshooting               | 24 |

## V

|                |    |
|----------------|----|
| Viewing height | 19 |
|----------------|----|

## W

|                  |    |
|------------------|----|
| Warning          |    |
| labels           | 10 |
| lights           | 10 |
| Warning labels   | 11 |
| Weight and Sizes | 28 |

## Z

|                          |    |
|--------------------------|----|
| ZEISS                    |    |
| Portal                   | 7  |
| Service agreements       | 22 |
| Zoom factor of objective | 15 |

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